

EASY-nLC Virtual Instrument for Xcalibur 2.1.x Quick Reference Guide

This guide describes how to install, configure, and use the Thermo EASY-nLC™ Virtual Instrument (VI) software with the Thermo Xcalibur™ data system version 2.1.x and the Thermo Foundation™ application version 1.0.2 or later. For information about Xcalibur specific operations, refer to the Xcalibur data system Help.

Contents

- Prerequisites
- Installation
- Removing a Previous Version
- Configuration
- Excluding the EASY-nLC from the VirusScan Enterprise
- Creating Methods in Xcalibur
- Creating Sequence Runs in Xcalibur
- Running a Sample Sequence
- Monitoring the Status
- Sample Information
- Additional Support
- Trademarks



Note When creating EASY-nLC methods and executing sample sequence runs, you must use the EASY-nLC Virtual Instrument (VI) application with the Xcalibur data system. For all other operations, such as configuration and maintenance tasks, use the touch monitor on the EASY-nLC instrument itself.

Prerequisites

Before installing the EASY-nLC Virtual Instrument software, ensure the following:

- Installation of Foundation version 1.0.2 or later
- Installation of Xcalibur version 2.1.x or later
- Administrator rights to install the EASY-nLC VI software for the Xcalibur application
- Removal of any earlier version of the EASY-nLC VI from the data system computer prior to installing the new version
(For guidance on how to remove the EASY-nLC VI software, see “[Removing a Previous Version](#)” on [page 2](#).)
- A network connection between the EASY-nLC instrument and the computer running the Xcalibur application
(To set up this connection, contact local IT support and consult the *EASY-nLC II User Guide*.)

IMPORTANT Avoid IP conflicts with other units on the network, especially the mass spectrometer. Refer to your mass spectrometer vendor for assistance in setting up the network connection between the host computer and additional peripheral units to the mass spectrometer.

- Access through port 6666 for connection to the EASY-nLC instrument.
- Connecting the EASY-nLC instrument and the mass spectrometer with the contact closure cable, and configuring the EASY-nLC system for one-way contact closure
(For further assistance, refer to the EASY-nLC or mass spectrometer manual as appropriate.)


Thermo Fisher Scientific is not responsible for problems that occur as a result of failure to comply with these instructions.

Installation

❖ To install the EASY-nLC Virtual Instrument software

1. Close all running programs before starting the installation. Remember to remove earlier versions of the EASY-nLC VI for the Xcalibur application before installing the new version of the software.
2. Load the LC Devices DVD.
3. Follow the instructions in the InstallShield Wizard.
4. On the Select Features page of the wizard, select the **Thermo** check box on the left, and then select the **Thermo EASY-nLC Xcalibur** check box on the right.
5. When the installation finishes, restart the data system computer.


The installation is complete when you log back in.

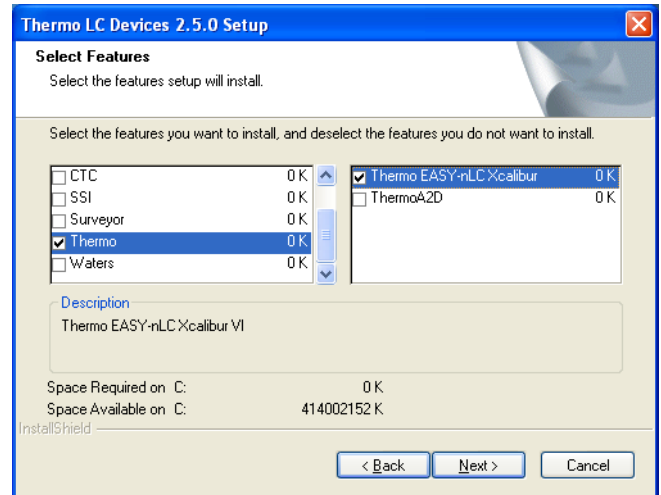

Removing a Previous Version

❖ To remove a previously installed version

1. Start **Add or Remove Programs** by choosing **Start > Settings > Control Panel**.
2. In the list of installed programs, locate **Thermo EASY-nLC Xcalibur VI**. Select it and click **Remove**.

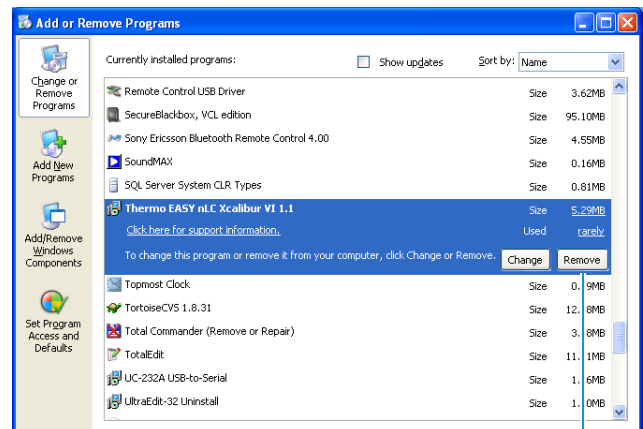

Remove button

Configuration

❖ To configure the EASY-nLC VI software for the Xcalibur data system

1. From your data system computer, choose **Start > Programs > Thermo Foundation 1.0 > Instrument Configuration**.

Note You can access Help at any time by pressing F1.

All available devices are listed in the Available Devices box; all configured devices are listed in the Configured Devices box.

2. Make sure that Device Types is set to **All** or **AS** (AutoSampler). Thermo EASY-nLC VI now appears in the Available Devices box.
The EASY-nLC VI for Xcalibur device does not appear if the Device Type is set to LC.
3. Select the **Thermo EASY-nLC** icon in the Available Devices box and click **Add** to move it to the Configured Devices box.

4. Select the **Thermo EASY-nLC** icon in the Configured Devices box and click **Configure**.

The Thermo EASY-nLC Configuration dialog box appears.

5. Type the IP address of the EASY-nLC system and your user name in the appropriate boxes.
You must have an account on the EASY-nLC system with the same user name.

The default user name is “admin,” but you can change this to any existing user account on the EASY-nLC system.

Note The EASY-nLC IP address is listed in the About box, which you can open by pressing the THERMO logo in the upper left corner of the touch screen.

The current user is listed in the lower left corner of the touch screen.

6. Click **Test Connection** to test the connection to the EASY-nLC system.

If the connection is established, the status box to the right of the Test Connection button changes to green (see the next figure). If a connection is not established, the status bar changes to red.

If an antivirus program or firewall is blocking the connection, the application displays this error message:



An established connection was aborted by the software in your host machine.

Contact your local IT administrator for assistance.

For information about excluding the EASY-nLC communication port from the McAfee virus scanner, see [“Excluding the EASY-nLC from the VirusScan Enterprise”](#) on page 5.


7. In the Autosampler Configuration area, select the autosampler model (**Autosampler ASA** or **Autosampler ASC**) and the plate type (**96-** or **384-well plates**, or a **4×6 (ASA)** or **6×8 (ASC)** vial holder).

Note All plate formats for the Autosampler ASC have six extra vial positions (1-V1 to 1-V6) on the side.

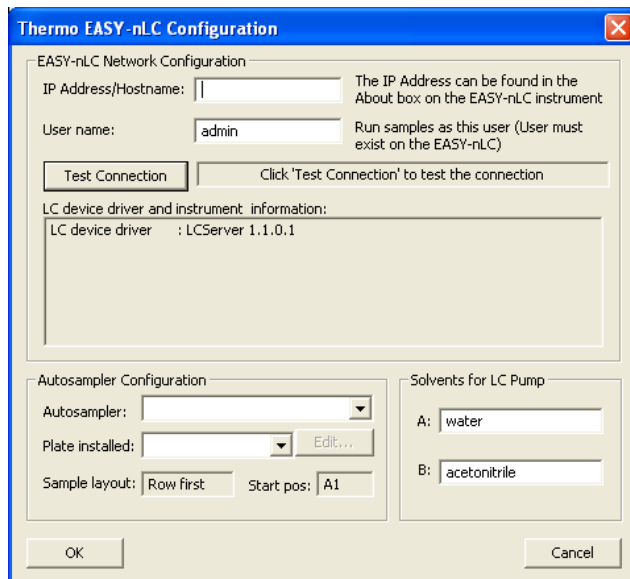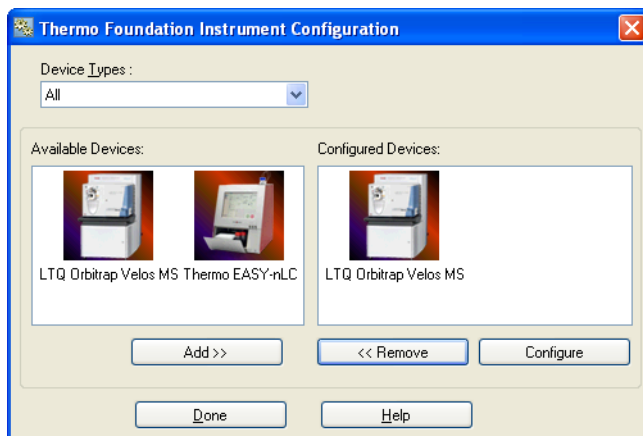

- In the Solvents for LC Pump area, type the names of the solvents in the boxes for solvent bottles A and B. Solvent bottle A typically contains LC/MS-grade water with 0.1% formic acid. Solvent bottle B typically contains LC/MS-grade acetonitrile with 0.1% formic acid.
- Configure additional parameters by clicking **Edit**.

The Thermo EASY-nLC Sample Position Selector dialog box opens.

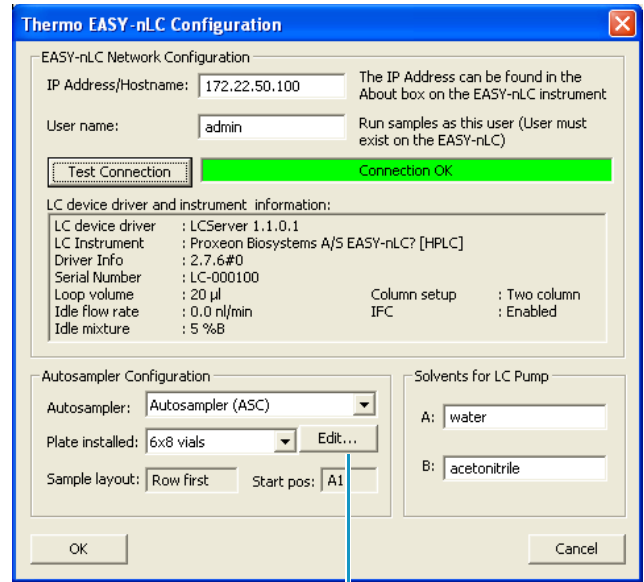


Edit button

- Select the default sample position (first sample to be run) by clicking a position in the Plate Rack area.

You can subsequently define a new default start position by clicking Clear All and then selecting a new position.

You can define the sample selection layout by selecting the Row First or Column First option in the lower left portion of the dialog box.


- Click **OK** to save your changes and exit the Thermo EASY-nLC Sample Position Selector dialog box.
- Click **OK** to exit the Thermo EASY-nLC Configuration dialog box.

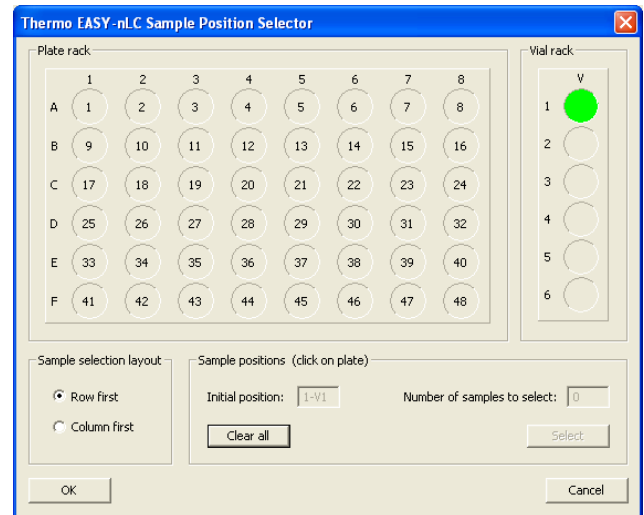

The McAfee™ virus scanner blocks the communication port for the Thermo EASY-nLC instrument when the data system computer uses the MicroSoft™ Windows™ 7 operating system.

Note For the Windows 7 operating system, you must exclude the EASY-nLC VI software from the McAfee blocking list. If your data system uses the Windows XP operating system, exclude the EASY-nLC VI software from the McAfee blocking list if your system is experiencing communication problems.

❖ **To exclude the EASY-nLC communication port from the VirusScan Enterprise 8.7.0i**

1. Log in to the data system computer as a user with administrative privileges.
2. Open the Virus Scan Console as follows:
 - a. Do one of the following:
 - On the Windows 7 taskbar, right-click the **McAfee** icon, .
 - or–
 - If the McAfee icon is hidden, click the Up arrow (on the left side of the notification area of the taskbar) to display the hidden icons. Then right-click the **McAfee** icon.

The McAfee shortcut menu appears.

- b. Choose **VirusScan Console**.

The VirusScan Console appears.

3. Right-click **Access Protection** and choose **Properties** from the shortcut menu.

The Access Protection Properties dialog box appears.

4. To open the Rules dialog box, do the following:
 - a. Under Categories, choose **Anti-virus Standard Protection**.
 - b. Under Rules, select **Prevent IRC Communication**.
 - c. Click **Edit**.

The Rules Details dialog box appears.

5. Modify these settings:
 - In the Processes to Include box, delete the asterisk.

Note When the Processes to Include box contains only a single asterisk, the virus scanner prevents IRC communication for all processes.

- In the Processes to Exclude box, type **Thermo***.

Note When the Processes to Exclude box contains the text Thermo*, the Access Protection service for Internet Relay Chat (IRC) communication processes ignores processes that begin with Thermo.

6. Click **OK** to close the Rules Details dialog box.
7. Click **OK** to close the Access Protection Properties dialog box and accept the new settings.
8. Close the VirusScan Console.

❖ **To create EASY-nLC methods in the Xcalibur data system**

1. Choose **GoTo > Instrument Setup**, or click the large **Instrument Setup** icon on the Xcalibur Roadmap Home Page. All installed devices are represented by icons in the left pane of the Instrument Setup window. The icon of your mass spectrometer will appear here.


Note

- Instrument methods (.meth file type) contain the chromatographic and data acquisition parameters for all the configured devices in your LC/MS system. Remember to define the parameters for all devices when creating new methods.
- The injection volume (sample pickup volume) is set in the method file, as it has to be coordinated with the rest of the method parameters, primarily the loading volume. You can override the injection volume in the Xcalibur sequence table; however, Thermo Fisher Scientific generally discourages doing so, having carefully considered all the method parameters.

2. Click the **Thermo EASY-nLC** icon and create your method.

The Xcalibur method editor is essentially similar to the method editor installed in the EASY-nLC software. For instructions on how to set up methods, refer to the *EASY-nLC II User Guide*.

3. Save the method before exiting the Instrument Setup window.

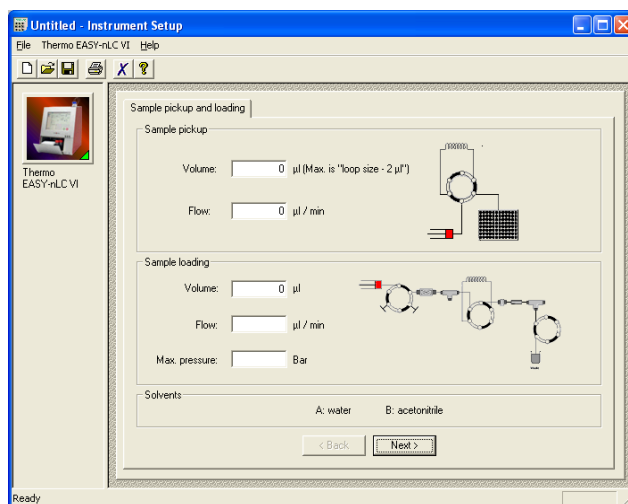


❖ To create sample sequence runs in the Xcalibur data system

1. Choose **View > Sequence Setup View** or click the large **Sequence Setup** icon on the Roadmap Home Page.
2. Choose **File > New**, or click the **New** icon in the toolbar to create a new sequence template.


For guidance on how to set up parameters in the New Sequence Template dialog box, refer to the Xcalibur user manual or press F1 to access the Help.

3. In the Samples area, click **Select Vials** to change the default well or vial selection.

The Thermo EASY-nLC Sample Position Selector dialog box appears.

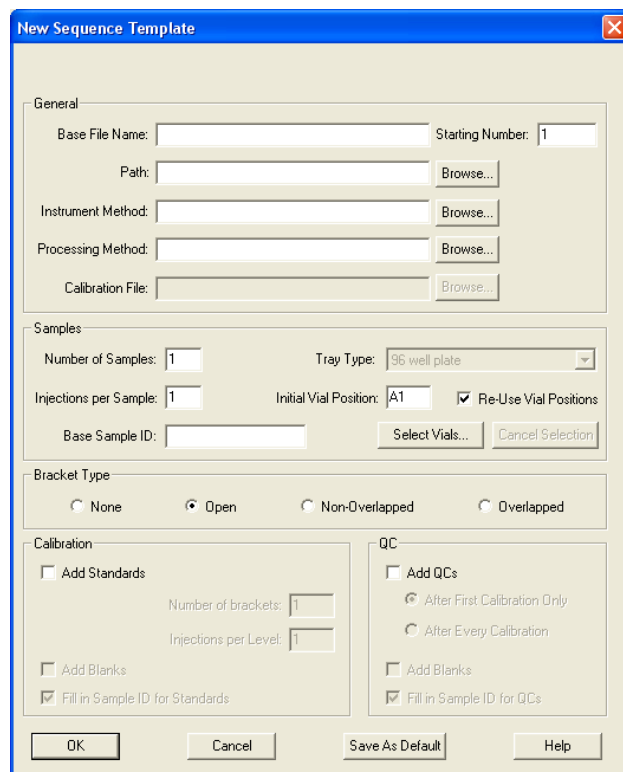

- Use the default start position, or define a new start position by clicking **Clear All** and then selecting a new position by clicking the plate.
- Use the default layout, can change the layout by selecting the **Row First** or **Column First** option.
- Type the number of additional samples to select in the Number of Samples to Select box and click **Select**. The consecutive wells and vials are then selected according to the option set in the Sample Selection Layout area.
- Click **OK** to return to the New Sequence Template dialog box.

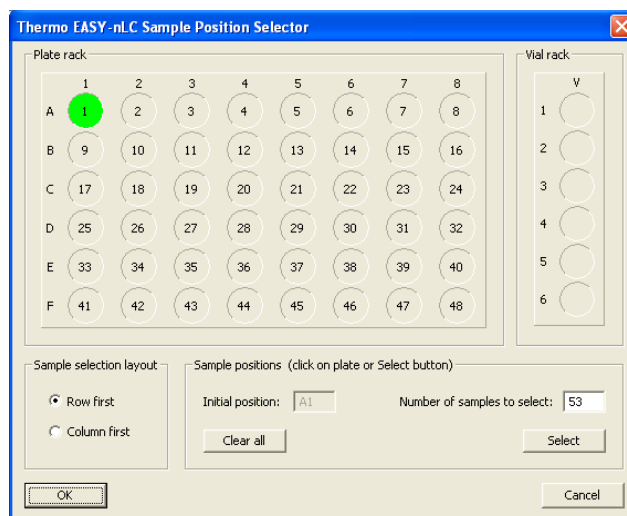

Note The position codes for the six extra vials on the side (Autosampler ASC) are 1-V1 to 1-V6.

- Click **OK** to create the sample sequence. All the sequence information is then listed in the Sequence Setup view on the Home Page.

	File Name	Path	Inst Meth	Position	Inj Vol
1	Assay01	C:\Xcalibur\Data\	C:\Xcalibur\methods\Protein Assay	A1	-1.0
2	Assay02	C:\Xcalibur\Data\	C:\Xcalibur\methods\Protein Assay	A2	-1.0
3	Assay03	C:\Xcalibur\Data\	C:\Xcalibur\methods\Protein Assay	A3	-1.0
4	Assay04	C:\Xcalibur\Data\	C:\Xcalibur\methods\Protein Assay	A4	-1.0
5	Assay05	C:\Xcalibur\Data\	C:\Xcalibur\methods\Protein Assay	B1	-1.0
6	Assay06	C:\Xcalibur\Data\	C:\Xcalibur\methods\Protein Assay	B2	-1.0
7	Assay07	C:\Xcalibur\Data\	C:\Xcalibur\methods\Protein Assay	B3	-1.0
8	Assay08	C:\Xcalibur\Data\	C:\Xcalibur\methods\Protein Assay	B4	-1.0
*					0.0

Note The sequence table contains a column called Inj Vol (injection volume). If the column does not appear, choose **Change > Column Arrangement** and add the Inj Vol column to the displayed columns. The value specified in this column overrules the value specified in the EASY-nLC method, as long as the value is between 0 and 48 μ L. This allows you to edit the injection volume of specific samples without changing the EASY-nLC method.

To ensure that the injection volume specified in the method file is used as intended, the default value in the Inj Vol column is set to -1 , an invalid injection volume. When the Xcalibur application detects the invalid value during a run, it automatically takes the actual injection volume from the corresponding method file, overwriting the -1 with the correct value.

If you choose to specify the injection volume manually in the Inj Vol column, ensure that the injection volume is sensible and complies with the size of the installed sample loop and other parameters in the method file (for example, load volume).

Never inject more than the *sample loop size* minus 2 μ L.

Running a Sample Sequence


❖ To run a single sample in the Xcalibur data system

1. Mark the specific row in the sequence table.
2. Choose **Actions > Run This Sample**, or click the **Run Sample** icon in the toolbar.

❖ To run a sequence in the Xcalibur data system

Choose **Actions > Run Sequence**, or click the **Run Sequence** icon in the toolbar.


Prior to execution of the sample or sequence run, the Run Sequence dialog box opens. For guidance on how to set up this dialog box, refer to the Xcalibur user manual or press F1 to access the Help.

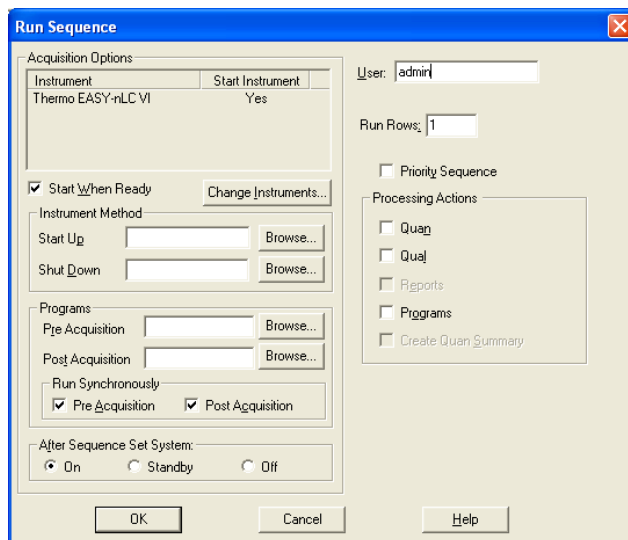

Monitoring the Status

❖ To monitor EASY-nLC status in the Xcalibur data system

- After installing and configuring the EASY-nLC VI software, and connecting the instrument to the Xcalibur data system, check the status in the Status pane of the Xcalibur Roadmap Home Page. A status message also appears.
- To check EASY-nLC method progress information and other EASY-nLC hardware readouts in real time, click **Ready To Download** at the top of the Status pane.

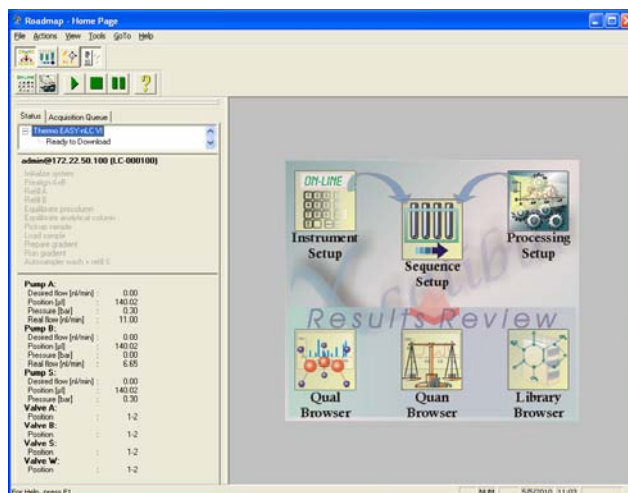

Sample Information

Using the Qual Browser application, you can access sample information and EASY-nLC method information in the raw MS data. Use any of these methods:

- On the Xcalibur Roadmap Home Page, click the **Qual Browser** icon.
- Open the MS raw file that you want to access method information from.
- Choose **View > Report > Sample Information** to view sample information.
- Choose **View > Report > Instrument Method** to view method information from all devices used to acquire the data (typically the EASY-nLC instrument and a mass spectrometer).
- Navigate through the pages using the arrow buttons in the toolbar or the arrow keys on the keyboard.

Additional Support

Visit www.thermofisher.com to find your local Thermo Fisher Scientific representative.

Trademarks

EASY-nLC and Foundation are trademarks and Xcalibur is a registered trademark of Thermo Fisher Scientific Inc. in the United States.

McAfee is a registered trademark of McAfee, Inc. in the United States and other countries.