

Investigation 4: Diffusion and Osmosis


Notes From the teacher

Day 1:

Before class: Read Learning Objectives through Procedure 1 and complete Day 1 Pre Lab.

Pre-Lab:

1. What is diffusion?
2. What is kinetic energy?
3. What environmental factors affect kinetic energy and diffusion?
4. Why are gradients important in diffusion?
5. Why are most cells small?
6. What is Brothelmal Blue?
7. Calculate the surface area of the following cube.
8. Calculate the volume of the following cube.

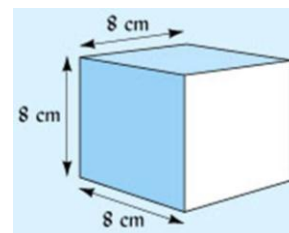

In class:

- Procedure 1, Step 2
In your lab notebook record sketches of your cubes. Label their dimensions. Provide the surface area, volume and ratio of each cube. Rank your 3 blocks fastest to slowest based on the speed at which you **think** the solution would diffuse through the entire block if it was placed in the solution.
- Procedure 1, DCYO
In your lab notebook follow the DCYO worksheet steps 2-4. *You do not have to write hypotheses or data tables.* Have your plan approved by Mrs. Enama.

Record your results as sketches of your cubes. Label their dimensions. Provide the surface area, volume, ratio and extent of diffusion of each cube.

Graph your results as the surface area to volume ratio versus the extent of diffusion.

Complete the analysis question in your lab notebook.

1. The size of some human cells is 0.01 mm. Using the formulas in this activity, calculate the surface to volume ratio of such a cell (assume 0.01 mm cube). Extrapolate the extent of diffusion into this living cell as compared to the smallest agar cube on your graph.

↻ Surface Area =
Length x width x # of sides

↻ Volume =
Length x width x height

↻ Surface Area Volume Ratio =
 $\frac{\text{Surface Area}}{\text{Volume}}$

↻ Extent of Diffusion =
 $\frac{\text{Total Cube Volume} - \text{Volume of cube that has not changed color}}{\text{Total Cube Volume}} \times 100$

Day 2

Before class: Read Procedure 2 and complete Day 2 Pre Lab.

Pre-Lab:

1. Why is it important for an IV solution to have salts in it?
2. What would happen if you were given pure water in your IV?
3. Describe how dialysis tubing is similar to a cell membrane.
4. Do you think dialysis tubing is permeable:
 - a. Water
 - b. Sucrose
 - c. NaCl
 - d. Glucose
 - e. Protein
5. Why is it important to leave room in your tubing when you knot it?
6. Calculate the percent change for the sample data provided.

Initial weight = 12.7g

Final weight = 13.6g

In class:

- Procedure 2, Steps 1-4
In your lab notebook record your set-up for each beaker (1-5). Make sure to include what solution is in the beaker and what solution is in the tubing.

As you perform the experiment record your results. (Results should include: initial weight, final weight, and percent change) For each beaker explain which way water moved, which way the solute(s) moved and why you think this occurred.

Complete the analysis question in your lab notebook.

1. From your results, which solutes, if any, diffused across the membrane and which, if any, were restricted? Why do you think this occurred?

Day 3:

Before Class: Read the Background Section, including water potential, Read Procedure 3, and complete Day 3 Pre Lab.

Pre-Lab

1. Describe the relative solute concentration and water potential in the following solutions:
 - a. Hypertonic
 - b. Hypotonic
2. How does water move in terms of:
 - a. Solute concentration
 - b. Water potential
3. Why do we need to consider pressure in plant cell osmosis but not animal cell osmosis?
4. What is the water potential in an open beaker of pure water?
5. Why does osmosis eventually stop in a plant cell, even if the environment is hypotonic?
6. Calculate the solute potential of a 0.1M NaCl solution at 25 C.
7. Calculate the solute potential of a 0.15M NaCl cell at 25 C.
8. Which way will the water move if the cell (#7) is placed in the solution (#6).
9. What must the turgor pressure equal if there is no net movement when the cell (#7) is placed in the solution (#6)?

In class:

- Review the lab protocol and water potential with the teacher. In your lab notebook follow the DCYO worksheet, do not write hypotheses. REMEMBER: the potatoes in solution will have to soak for 24 hours. Have your plan approved by the teacher.

Day 4/5

Before class: N/A

In class:

- Conduct your investigation
In your lab notebook record your results. (Results should include: color of solution, potato's initial weight, potato's final weight, potato's percent change, and which way the water moved.

Complete the analysis question in your lab notebook.

1. Which color solution had the highest concentration of sucrose? How do you know this?
2. Match each color solution to its actual sucrose concentration. The concentrations were 0.2 M, 0.4 M, 0.6 M, 0.8 M, and 1.0 M.

3. Once the color solutions have been matched to their actual sucrose concentrations calculate the water potential of the potato core in each solution. Show your calculations and explain your reasoning.
4. If you looked at your potato cores under the microscope, describe what you think you would see in each color solution.

Cellular Processes: Energy and Communication

INVESTIGATION 4

DIFFUSION AND OSMOSIS

What causes my plants to wilt if I forget to water them?


■ BACKGROUND

Cells must move materials through membranes and throughout cytoplasm in order to maintain homeostasis. The movement is regulated because cellular membranes, including the plasma and organelle membranes, are selectively permeable. Membranes are phospholipid bilayers containing embedded proteins; the phospholipid fatty acids limit the movement of water because of their hydrophobic characteristics.

The cellular environment is aqueous, meaning that the solvent in which the solutes, such as salts and organic molecules, dissolve is water. Water may pass slowly through the membrane by osmosis or through specialized protein channels called aquaporins. Aquaporins allow the water to move more quickly than it would through osmosis. Most other substances, such as ions, move through protein channels, while larger molecules, including carbohydrates, move through transport proteins.

The simplest form of movement is diffusion, in which solutes move from an area of high concentration to an area of low concentration; diffusion is directly related to molecular kinetic energy. Diffusion does not require energy input by cells. The movement of a solute from an area of low concentration to an area of high concentration requires energy input in the form of ATP and protein carriers called pumps.

Water moves through membranes by diffusion; the movement of water through membranes is called osmosis. Like solutes, water moves down its concentration gradient. Water moves from areas of high potential (high free water concentration) and low solute concentration to areas of low potential (low free water concentration) and high solute concentration. Solute molecules decrease the concentration of free water, since water molecules surround the solute molecules. The terms *hypertonic*, *hypotonic*, and *isotonic* are used to describe solutions separated by selectively permeable membranes. A hypertonic solution has a higher solute concentration and a lower water potential as compared to the other solution; therefore, water will move into the hypertonic solution through the membrane by osmosis. A hypotonic solution has a lower solute concentration and a higher water potential than the solution on the other side of the membrane; water will move down its concentration gradient into the other solution. Isotonic solutions have equal water potentials.



In nonwalled cells, such as animal cells, the movement of water into and out of a cell is affected by the relative solute concentration on either side of the plasma membrane. As water moves out of the cell, the cell shrinks; if water moves into the cell, it swells and may eventually burst. In walled cells, including fungal and plant cells, osmosis is affected not only by the solute concentration, but also by the resistance to water movement in the cell by the cell wall. This resistance is called turgor pressure. The presence of a cell wall prevents the cells from bursting as water enters; however, pressure builds up inside the cell and affects the rate of osmosis.

Water movement in plants is important in water transport from the roots into the shoots and leaves. You likely will explore this specialized movement called transpiration in another lab investigation.

■ Understanding Water Potential

Water potential predicts which way water diffuses through plant tissues and is abbreviated by the Greek letter psi (ψ). Water potential is the free energy per mole of water and is calculated from two major components: (1) the solute potential (ψ_s), which is dependent on solute concentration, and (2) the pressure potential (ψ_p), which results from the exertion of pressure—either positive or negative (tension) — on a solution. The solute potential is also called the osmotic potential.

$$\psi = \psi_p + \psi_s$$

Water Potential = Pressure Potential + Solute Potential

Water moves from an area of higher water potential or higher free energy to an area of lower water potential or lower free energy. Water potential measures the tendency of water to diffuse from one compartment to another compartment.

The water potential of pure water in an open beaker is zero ($\psi = 0$) because both the solute and pressure potentials are zero ($\psi_s = 0$; $\psi_p = 0$). An increase in positive pressure raises the pressure potential and the water potential. The addition of solute to the water lowers the solute potential and therefore decreases the water potential. This means that a solution at atmospheric pressure has a negative water potential due to the solute.

The solute potential (ψ_s) = $-iCRT$, where i is the ionization constant, C is the molar concentration, R is the pressure constant ($R = 0.0831$ liter bars/mole-K), and T is the temperature in K ($273 + ^\circ\text{C}$).

A 0.15 M solution of sucrose at atmospheric pressure ($\psi_p = 0$) and 25°C has an osmotic potential of -3.7 bars and a water potential of -3.7 bars. A bar is a metric measure of pressure and is the same as 1 atmosphere at sea level. A 0.15 M NaCl solution contains 2 ions, Na^+ and Cl^- ; therefore $i = 2$ and the water potential = -7.4 bars.

When a cell's cytoplasm is separated from pure water by a selectively permeable membrane, water moves from the surrounding area, where the water potential is higher ($\psi = 0$), into the cell, where water potential is lower because of solutes in the cytoplasm

(ψ is negative). It is assumed that the solute is not diffusing (Figure 1a). The movement of water into the cell causes the cell to swell, and the cell membrane pushes against the cell wall to produce an increase in pressure. This pressure, which counteracts the diffusion of water into the cell, is called turgor pressure.

Over time, enough positive turgor pressure builds up to oppose the more negative solute potential of the cell. Eventually, the water potential of the cell equals the water potential of the pure water outside the cell (ψ of cell = ψ of pure water = 0). At this point, a dynamic equilibrium is reached and net water movement ceases (Figure 1b).



Figures 1a-b. Plant cell in pure water. The water potential was calculated at the beginning of the experiment (a) and after water movement reached dynamic equilibrium and the net water movement was zero (b).

If solute is added to the water surrounding the plant cell, the water potential of the solution surrounding the cell decreases. If enough solute is added, the water potential outside the cell is equal to the water potential inside the cell, and there will be no net movement of water. However, the solute concentrations inside and outside the cell are not equal, because the water potential inside the cell results from the combination of both the turgor pressure (ψ_p) and the solute pressure (ψ_s). (See Figure 2.)

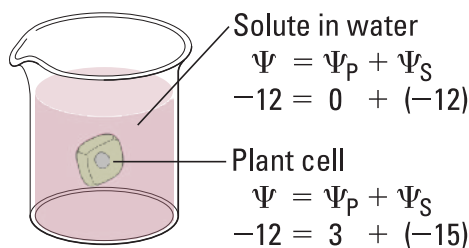



Figure 2. Plant cell in an aqueous solution. The water potential of the cell equals that of surrounding solution at dynamic equilibrium. The cell's water potential equals the sum of the turgor pressure potential plus the solute potential. The solute potentials of the solution and of the cell are not equal.

If more solute is added to the water surrounding the cell, water will leave the cell, moving from an area of higher water potential to an area of lower water potential. The water loss causes the cell to lose turgor. A continued loss of water will cause the cell membrane to shrink away from the cell wall, and the cell will plasmolyze.

- 
- Calculate the solute potential of a 0.1 M NaCl solution at 25°C. If the concentration of NaCl inside the plant cell is 0.15 M, which way will the water diffuse if the cell is placed into the 0.1 M NaCl solutions?
 - What must the turgor pressure equal if there is no net diffusion between the solution and the cell?

■ Learning Objectives

- To investigate the relationship among surface area, volume, and the rate of diffusion
- To design experiments to measure the rate of osmosis in a model system
- To investigate osmosis in plant cells
- To design an experiment to measure water potential in plant cells
- To analyze the data collected in the experiments and make predictions about molecular movement through cellular membranes
- To work collaboratively to design experiments and analyze results
- To connect the concepts of diffusion and osmosis to the cell structure and function

■ General Safety Precautions

You must wear safety glasses or goggles, aprons, and gloves because you will be working with acids and caustic chemicals. The HCl and NaOH solutions will cause chemical burns, and you should use these solutions in spill-proof trays or pans. Follow your teacher's instructions carefully. Do not work in the laboratory without your teacher's supervision. Talk to your teacher if you have any questions or concerns about the experiments.

■ THE INVESTIGATIONS

This investigation consists of three parts. In Procedure 1, you use artificial cells to study the relationship of surface area and volume. In Procedure 2, you create models of living cells to explore osmosis and diffusion. You finish by observing osmosis in living cells (Procedure 3). All three sections of the investigation provide opportunities for you to design and conduct your own experiments.

■ Getting Started

These questions are designed to help you understand kinetic energy, osmosis, and diffusion and to prepare for your investigations.

- What is kinetic energy, and how does it differ from potential energy?
- What environmental factors affect kinetic energy and diffusion?

- How do these factors alter diffusion rates?
- Why are gradients important in diffusion and osmosis?
- What is the explanation for the fact that most cells are small and have cell membranes with many convolutions?
- Will water move into or out of a plant cell if the cell has a higher water potential than the surrounding environment?
- What would happen if you applied saltwater to a plant?
- How does a plant cell control its internal (turgor) pressure?

■ Procedure 1: Surface Area and Cell Size

Cell size and shape are important factors in determining the rate of diffusion. Think about cells with specialized functions, such as the epithelial cells that line the small intestine or plant root hairs.

- What is the shape of these cells?
- What size are the cells?
- How do small intestinal epithelial and root hair cells function in nutrient procurement?

Materials


- | | |
|---|---|
| • 2% agar containing NaOH and the pH indicator dye phenolphthalein | disposable plates); unserrated knives; or scalpels from dissection kits |
| • Vinegar Solution | • Metric rulers |
| • 0.1 M HCl | • Petri dishes and test tubes |
| • 0.1 M NaOH | • Brothemal Blue agar preparation |
| • Squares of hard, thin plastic (from | |

Step 1 ~~Place some phenolphthalein in two test tubes. Add 0.1 M HCl to one test tube, swirl to mix the solutions, and observe the color. Using the same procedure, add 0.1 M NaOH to the other test tube. Remember to record your observations as you were instructed.~~

- ~~Which solution is an acid?~~
- ~~Which solution is a base?~~
- ~~What color is the dye in the base? In the acid?~~
- ~~What color is the dye when mixed with the base?~~

Skip Part 1 - We will be using Brothemal Blue Agar. Brothemal Blue is an acid indicator.

It is blue and turns yellow in the presence of acid.



Step 2 Using a dull knife or a thin strip of hard plastic, cut three blocks of agar of different sizes. Please keep your agar cubes smaller than 2cm per side.

These three blocks will be your models for cells.

- What is the surface area of each of your three cells?
- What is the total volume of each of your cells?
- If you put each of the blocks into a solution, into which block would that solution diffuse throughout the entire block fastest? Slowest? How do you explain the difference?

Alternative Method

~~Mix one packet of unflavored gelatin with 237 mL of water; add 2.5 mL 1% phenolphthalein and a few drops of 0.1 M NaOH. The solution should be bright pink. Pour the gelatin mixture into shallow pans and refrigerate overnight.~~

~~You may use white vinegar in place of the 0.1 M HCl.~~

Designing and Conducting Your Investigation

Using the materials listed earlier, design an experiment to test the predictions you just made regarding the relationship of surface area and volume in the artificial cells to the diffusion rate using the Brothelmal Blue Agar and Vinegar solution. NOTE: you will need your blocks to soak for 20 to 30 minutes. Once you have finished planning your experiment, have your teacher check your design. When you have an approved design, run your experiment and record your results. Do your experimental results support your predictions?

Procedure 2: Modeling Diffusion and Osmosis

You are in the hospital and need intravenous fluids. You read the label on the IV bag, which lists all of the solutes in the water.

- Why is it important for an IV solution to have salts in it?
- What would happen if you were given pure water in an IV?
- How would you determine the best concentration of solutes to give a patient in need of fluids *before* you introduced the fluids into the patient's body?

In this experiment, you will create models of living cells using dialysis tubing. Like cell membranes, dialysis tubing is made from a material that is selectively permeable to water and some solutes. You will fill your model cells with different solutions and determine the rate of diffusion.

Materials

- Distilled or tap water
 - 1 M sucrose
 - 1 M NaCl
 - 1 M glucose
 - 5% ovalbumin (egg white protein)
 - 20 cm-long dialysis tubing
 - Cups
 - Balances
- How can you use weights of the filled cell models to determine the rate and direction of diffusion? What would be an appropriate control for the procedure you just described?
 - Suppose you could test other things besides weights of the dialysis tubes. How could you determine the rates and directions of diffusion of water, sucrose, NaCl, glucose, and ovalbumin?
 - Will protein diffuse? Will it affect the rate of diffusion of other molecules?

Step 1 Choose up to four pairs of different solutions. One solution from each pair will be in the model cell of dialysis tubing, and the other will be outside the cell in the cup. Your fifth model cell will have water inside and outside; this is your control. Before starting, use your knowledge about solute gradients to predict whether the water will diffuse into or out of the cell. Make sure you label the cups to indicate what solution is inside the cell and inside the cup.

Step 2 Make dialysis tubing cells by tying a knot in one end of five pieces of dialysis tubing. Fill each “cell” with 10 mL of the solution you chose for the inside, and knot the other end, leaving enough space for water to diffuse into the cell.

Step 3 Weigh each cell, record the initial weight, and then place it into a cup filled with the second solution for that pair. Weigh the cell after 30 minutes and record the final weight.

Step 4 Calculate the percent change in weight using the following formula:
 $(\text{final} - \text{initial}) / \text{initial} \times 100$. Record your results.

- Which pair(s) that you tested did not have a change in weight? How can you explain this?
- If you compared 1 M solutions, was a 1 M NaCl solution more or less hypertonic than a 1 M sucrose solution? What is your evidence? What about 1 M NaCl and 1 M glucose and 1 M sucrose?
- Does the protein solution have a high molarity? What is evidence for your conclusion?
- How could you test for the diffusion of glucose?
- Based on what you learned from your experiment, how could you determine the solute concentration inside a living cell?

■ ~~Designing and Conducting Your Investigation~~



~~Living cell membranes are selectively permeable and contain protein channels that permit the passage of water and molecules. In some respects, the dialysis tubing you used is similar to a cell membrane, and you can use it to explore osmosis in greater depth. Think about the questions that came up as you worked through the investigation. What unanswered questions do you still have about osmosis that you could investigate further?~~

~~Using the available materials, design an investigation to answer one of your questions. Have your teacher check your design first. Remember to record your results, and be sure to use appropriate controls.~~

~~These questions can help jump start your thinking.~~

- ~~• What factors determine the rate and direction of osmosis?~~
- ~~• What would you predict if you used a starch solution instead of the protein?~~
- ~~• Can you diagram the flow of water based upon the contents of your model cell and the surrounding solution?~~
- ~~• When will the net osmosis rate equal zero in your model cells? Will it ever truly be zero?~~
- ~~• Based upon your observations, can you predict the direction of osmosis in living cells when the cells are placed in various solutions?~~
- ~~• How is the dialysis tubing functionally different from a cellular membrane?~~

■ Procedure 3: Observing Osmosis in Living Cells

The interactions between selectively permeable membranes, water, and solutes are important in cellular and organismal functions. For example, water and nutrients move from plant roots to the leaves and shoots because of differences in water potentials. Based upon what you know and what you have learned about osmosis, diffusion, and water potential in the course of your investigations, think about these questions.

- What would happen if you applied saltwater to the roots of a plant? Why?
- What are two different ways a plant could control turgor pressure, a name for internal water potential within its cells? Is this a sufficient definition for turgor pressure?
- Will water move into or out of a plant cell if the cell has a higher water potential than its surrounding environment?

Step 1 ~~Start by looking at a single leaf blade from either *Elodea* (a water plant) or a leaf-like structure from *Mnium hornum* (a moss) under the light microscope. If you need assistance, your teacher will show you how to place specimens on a slide.~~

- ~~• Where is the cell membrane in relation to the cell wall? Can you see the two structures easily? Why or why not?~~
- ~~• What parts of the cell that you see control the water concentration inside the cell?~~

~~Back in Procedure 2 you tested diffusion and osmosis properties of several solutions. Now you are going to determine how they affect plant cell turgor pressure.~~

- ~~• What changes do you expect to see when the cells are exposed to the solutions?~~
- ~~• How will you know if a particular treatment is increasing turgor pressure? If it is reducing turgor pressure?~~
- ~~• How could you determine which solution is isotonic to the cells?~~

Step 2

Click the following link if you do not remember what a plant cell looks like

undergoing plasmolysis: <http://www.dnatube.com/video/2775/Onion-Cells-Plasmolysis>

■ Designing and Conducting Your Investigation

- Potatoes, sweet potatoes, or yams
- Cork borers or french fry cutter
- Balances
- Metric rulers
- Cups
- Color-coded sucrose solutions of different, but unlabeled, concentrations prepared by your teacher

Design an experiment to identify the concentrations of the sucrose solutions and use the solutions to determine the water potential of the plant tissues. (You might want to review the information on water potential described in Understanding Water Potential.)

Use the following questions to guide your investigation:

- How can you measure the plant pieces to determine the rate of osmosis?
- How would you calculate the water potential in the cells?
- Which solution had a water potential equal to that of the plant cells? How do you know?
- Was the water potential in the different plants the same?
- How does this compare to your previous determinations in the *Elodea* cells?
- What would your results be if the potato were placed in a dry area for several days before your experiment?
- When potatoes are in the ground, do they swell with water when it rains? If not, how do you explain that, and if so, what would be the advantage or disadvantage?



■ Analyzing Results

1. Why are most cells small, and why do they have cell membranes with many convolutions?
2. What organelles inside the cell have membranes with many convolutions? Why?
3. Do you think osmosis occurs when a cell is in an isotonic solution? Explain your reasoning.

■ Where Can You Go from Here?

Do you think that fungal cells have turgor pressure? Design an experiment to test your hypothesis.