



DNA MODIFYING ENZYMES

Gibson Assembly<sup>®</sup> Master Mix

Instruction Manual

NEB #E2611S/L  
10/50 reactions

 NEW ENGLAND  
**BioLabs**<sup>®</sup> Inc.  
*enabling technologies in the life sciences*



**Limited Use Label License:** This Product and its components and products therefrom, are to be used for internal research purposes for the sole benefit of the purchaser. They may not be used for any other purpose, including, but not limited to, use in drugs, diagnostics, therapeutics or in humans. This Product and its components and products therefrom may not be transferred or sold to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service of any kind to third parties, including, without limitation, reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact SGI-DNA at [licensing@sgidna.com](mailto:licensing@sgidna.com).

**Limited Warranty:** The Gibson Assembly Master Mix is warranted to perform according to the specifications stated on the certificate of analysis. No other warranty is made, whether express or implied, including any warranty of merchantability or fitness for a particular purpose. This warranty limits NEB's and its licensors' liability to only the price of the kit. Neither NEB nor its licensors shall have any responsibility or liability for any special, incidental, indirect or consequential loss or damage whatsoever.

**Limited Label License:** This product is covered by certain patents outside the United States (the "Non-US Patents") corresponding to the expired U.S. Patent No. 5,436,149 (claims 6-16). The purchase of this product conveys to the purchaser only the limited, non-transferable right under the Non-US Patents to use the purchased quantity of the product for the purchaser's own research by the purchaser only. No rights are granted to the purchaser with respect to the Non-US Patents to sell, modify for resale or otherwise transfer this product, either alone or as a component of another product, to any third party. Takara Bio reserves all other rights to the Non-US Patents, and this product may not be used in any manner other than as provided herein. For information on obtaining a license with respect to the Non-US Patents to use this product for purposes other than research, please contact Takara Bio Inc., Seto 3-4-1, Otsu, Shiga 520-2193, Japan (Fax +81-77-543-9254).


NEW ENGLAND BIOLABS® and Q5® are registered trademarks of New England Biolabs, Inc.

GIBSON ASSEMBLY® is a registered trademark of Synthetic Genomics, Inc.

MILLI-Q® is a registered trademark of Millipore, Inc.

NANODROP™ is a trademark of Thermo Fisher Scientific, Inc.

SYNTHETIC GENOMICS® is a registered trademark of Synthetic Genomics, Inc.

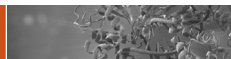

## Table of Contents:

---

Introduction.....	2
Specification.....	3
Overview of Gibson Assembly Master Mix Protocol.....	3
Design and PCR Amplification of Fragments for Gibson Assembly.....	3
Gibson Assembly Reaction.....	11
Chemically Competent Cells Transformation Protocol.....	12
Electrocompetent Cells Transformation Protocol.....	12
Usage Notes.....	13
Frequently Asked Questions (FAQs).....	14
Troubleshooting.....	17
References.....	19
Ordering Information.....	20

## Components:

---

*Store at -20°C. Thaw, vortex thoroughly before use and keep on ice.*

Gibson Assembly Master Mix (2X)

Positive Control

2 overlapping dsDNA fragments for control assembly.

### Required Materials Not Included:

DNA Polymerases (for generating PCR products):

We recommend Q5® High-Fidelity DNA Polymerase (NEB #M0491) or related products, such as Q5 Hot Start Flex DNA Polymerase (NEB #M0493), Q5 Hot Start Flex 2X Master Mix (NEB #M0494).

LB (Luria-Bertani) plates with appropriate antibiotic.

SOC Outgrowth Medium (NEB #B9020).

Competent Cells:

We recommend NEB 5-alpha Competent *E. coli* (High Efficiency, NEB #C2987).

For assembled products greater than 10 kb, NEB recommends using NEB 10-beta

Competent *E. coli* (High Efficiency, NEB #C3019) or NEB 10-beta Electrocompetent *E. coli* (NEB #C3020).

## Introduction:


Gibson Assembly was developed by Dr. Daniel Gibson and his colleagues at the J. Craig Venter Institute and licensed to NEB by Synthetic Genomics, Inc. It allows for successful assembly of multiple DNA fragments, regardless of fragment length or end compatibility. It has been rapidly adopted by the synthetic biology community due to its ease-of-use, flexibility and suitability for large DNA constructs.

Gibson Assembly efficiently joins multiple overlapping DNA fragments in a single-tube isothermal reaction (1,2). The Gibson Assembly Master Mix includes three different enzymatic activities that perform in a single buffer:

- The exonuclease creates single-stranded 3' overhangs that facilitate the annealing of fragments that share complementarity at one end (overlap region).
- The proprietary polymerase fills in gaps within each annealed fragment.
- The DNA ligase seals nicks in the assembled DNA.

The end result is a double-stranded fully sealed DNA molecule that can serve as template for PCR, RCA or a variety of other molecular biology applications, including direct transformation. The method has been successfully used by Gibson's group and others to assemble oligonucleotides, DNA with varied overlaps (15–80 bp) and fragments hundreds of kilobases long (1–2).

Figure 1: Overview of the Gibson Assembly Method

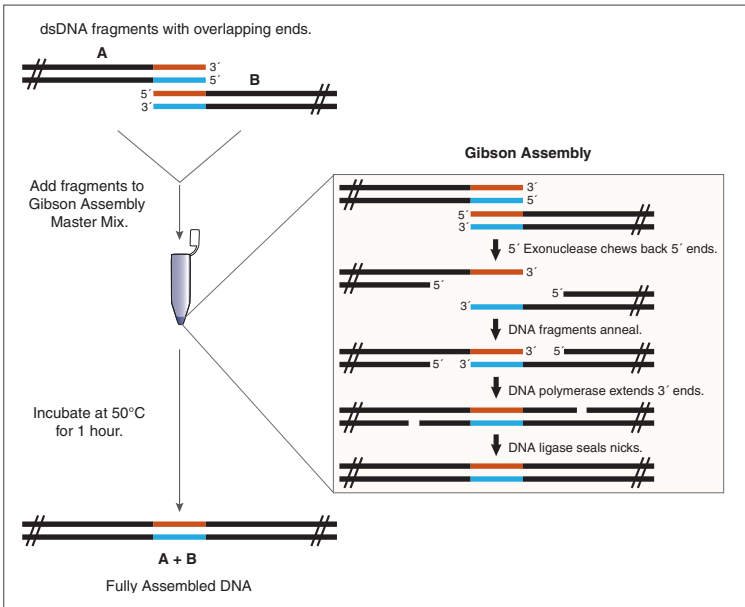

## Specification:

10  $\mu$ l of 2X Gibson Assembly Master Mix was incubated with 6 fragments (5 fragments of 400 bp and one of 2,780 bp, with 40 bp overlap, 0.05 pmol each) in a final volume of 20  $\mu$ l at 50°C for 60 minutes. NEB 5-alpha Competent *E. coli* (NEB #C2987) were transformed with 2  $\mu$ l of the master mix/fragment mixture using the transformation protocol on page 11. Greater than 100 white colonies were observed when 1/10 of the outgrowth was spread on an ampicillin plate with IPTG/Xgal and incubated overnight.

## Overview of Gibson Assembly Master Mix Protocol:

- Design primers to amplify fragments (and/or vector) with appropriate overlaps (see pages 3–7)
- PCR amplify fragments and/or vector using a high fidelity DNA polymerase (DNA can also be prepared using a restriction digest)
- Confirm and determine concentration of fragments using agarose gel electrophoresis, a Nanodrop™ instrument or other method
- Add DNAs to Gibson Assembly Master Mix and incubate at 50°C for 1 hour
- Transform into *E. coli* or use directly in other applications

## Design and PCR Amplification of Fragments for Gibson Assembly:

*Note: We highly recommend using our web tool, NEBuilder™, available at [www.NEBGibson.com](http://www.NEBGibson.com), to design PCR primers with overlapping sequences between the adjacent DNA fragments and for their assembly into a cloning vector.*



NEBuilder is the fastest and easiest approach to obtain ready-to-use sequences for overlapping primers. However, it does not give details about the primer design workflow. In some cases, it might be appropriate to further manually alter primer sequences in order to adapt them for the use in more complex assemblies, such as those that include site-specific mutagenesis. For this purpose, it is absolutely necessary to understand the general requirements and rules that apply to PCR primers used in conjunction with Gibson Assembly. The sections below offer step-by-step directions and recommendations for the manual design of primers for the assembly of two or more PCR fragments, as well as primer design for assembly of PCR fragments into a cloning vector prepared either by PCR or by restriction digestion.

## Structure of the Overlapping Primers

PCR primers for use in Gibson Assembly must have two sequence components:

- an overlap sequence, required for the assembly of adjacent fragments;
- a gene-specific sequence, required for template priming during PCR amplification;

The non-priming overlap sequence is added at the 5'-end of the primer. This sequence is homologous to the 5'-terminal sequence of the adjacent fragment. The length of overlap sequence is dependent on the GC content of the sequences.

The priming gene-specific sequence is added at the 3'-end of the primer after the overlap sequence. The priming sequence should meet the criteria required for template annealing during PCR amplification.



The  $T_m$  of the 3' gene-specific sequence of the primer can be calculated using the  $T_m$  calculator found on the NEB website at <http://www.neb.com/TmCalculator>.

## General Recommendations for Design of Overlapping Primers

To achieve efficient assembly of PCR fragments into a vector, we suggest using a 15–25 nt overlap with a  $T_m$  equal to or greater than 48°C (assuming A-T pair = 2°C and G-C pair = 4°C, Fig. 2A, Step I). To prevent errors in primer design it is highly recommended to first perform DNA fragment assembly *in silico* and create a final sequence file displaying both DNA strands. This virtual sequence may then be used as a template to design overlapping primers. Figure 2A shows the workflow for overlapping primer design by using an *in silico*-created DNA sequence file. First, mark the junctions between the adjacent fragments 1, 2 and 3 (Fig. 2A, step II). Next, at or near each junction choose 15–25 nucleotide sequences to serve as the overlap region between the two adjacent fragments (Fig. 2A, step III). For the best fit, in terms of length and  $T_m$ , the overlap sequence can be composed of nucleotides which belong to only one fragment (overlap shown in blue) or it can be split between the two adjacent fragments in any combination (overlap shown in orange). Mark the first 5' and the last 3' nucleotide of the overlap sequence on both DNA strands (boxed sequence). Finally, starting from the first 5' nucleotide, copy the entire overlap sequence in the 5' to 3' direction and, if necessary, continue to add additional nucleotides to the 3' end until the gene-specific priming sequence length is reached (Fig. 2A, step IV). The reverse overlapping primer is designed following the same steps as described above but copying the sequence from the complementary DNA strand in the 5' to 3' direction. Keep in mind that the two primers sharing the same overlap sequence are always used in separate PCR reactions, each in combination with the primer, which primes the complementary sequence on the opposite end of the respective DNA fragment (Fig. 2A, step V).

Figure 2A: Primer design using *in silico* created final DNA sequence file.

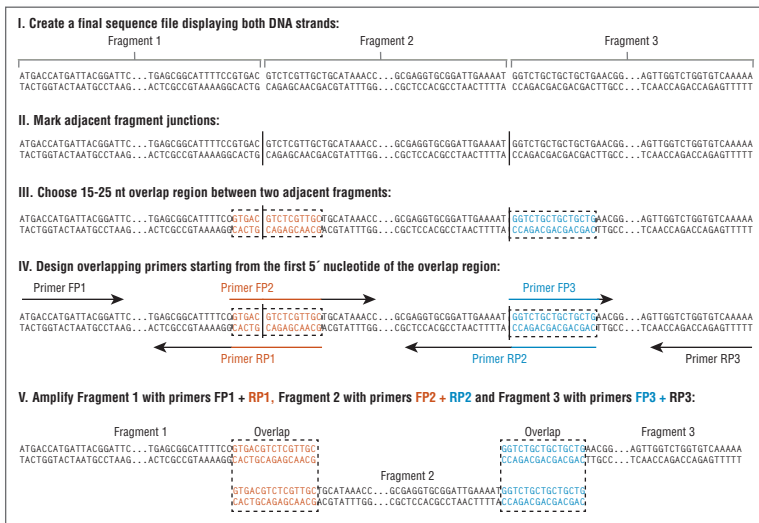


Figure 2B: Primer design for PCR-generated vector and insert using *in silico* created final DNA sequence file.



## Primer Design for PCR-Generated Vector and Insert


For the purposes of primer design, the vector and the insert may be viewed as two PCR fragments that have to be assembled into a circular DNA molecule. This means that the primer design rules described above may also be applied for generation of the vector fragment and the insert fragment sharing overlapping ends. Use the *in silico*-created final sequence file as a template to design overlapping primers between the vector and the insert by accomplishing the same steps as described above, as shown in Figure 2B.

If you intend to use PCR-generated vector for one specific insertion, then the overlap sequence may be split between the vector and the insert in any combination to make shorter primers (Figure 2B, step III, overlap shown in orange). However, if the same PCR-generated vector will be used for assembly of various inserts, then the entire overlap sequence must originate from the vector sequence and must be added to primers that will be used to amplify the insert (Figure 2B, step III, overlap shown in blue). The latter case is also illustrated in Figure 3 for assembly of the *lacZ* gene into a pET21a vector. The pET21a forward primer (orange arrow) and the reverse primer (blue arrow) start at the position where the *lacZ* gene must be inserted. Both vector-specific primers completely match the vector sequence on the respective strands marked in orange and blue. This inverse PCR strategy yields a linear vector fragment. Generally, 10–100 pg of a vector is recommended as a template in the inverse PCR reaction.

To amplify the *lacZ* gene, both forward and reverse *lacZ*-specific priming sequences (gray) at their 5' end are fused with 20-nt vector sequences to be used as overlap sequences in assembly with the vector. Within the *lacZ* Forward PCR primer, the overlap sequence (orange) is identical to the 20-nt terminal sequence on the top strand (orange) of the vector's left-arm (in the 5'→3' direction). Within the *lacZ* Reverse PCR primer, the overlap sequence (blue) is identical to the 21-nt terminal sequence on the bottom strand (blue) of the vector's right-arm (in the 5'→3' direction). The length of the overlap sequence is determined by the number of nucleotides needed to reach a  $T_m \geq 48^\circ\text{C}$ . If necessary, one may add additional nucleotides between the overlap sequence and the *lacZ*-specific sequence, for example, to introduce a unique restriction site.



Figure 3: Primer Design for Vector pET21a and *lacZ* Gene Assembly.

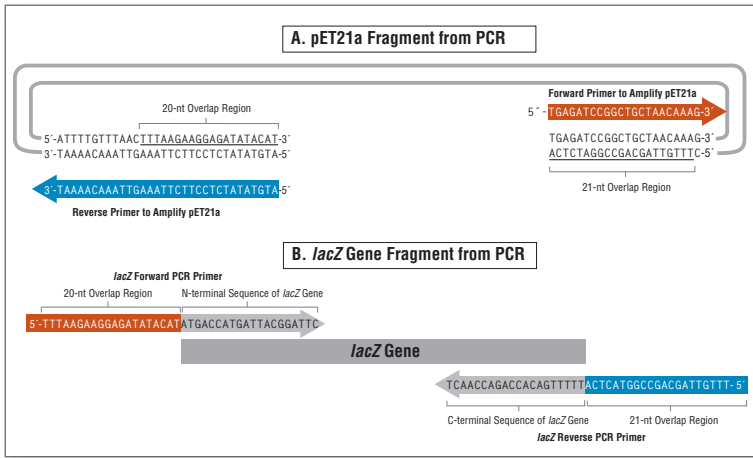

## Primer Design for Assembly of Restriction Enzyme Digested Vector and PCR-Generated Insert

Restriction enzyme-treated vectors can have 5'-overhangs, 3'-overhangs or blunt ends. When vector is linearized by restriction digestion, the entire overlap sequence must originate from the vector sequence and must be added to primers that will be used to amplify the insert. The overlap region of the forward primer for the gene of interest (orange) should line up with the 3' end of the overhang on the vector's left arm and extend back until the  $T_m \geq 48^\circ\text{C}$  (Fig.4A, Left side). This primer also includes gene-specific sequence at the 3'-end (gray). Keep in mind that the restriction site, which was used to digest the vector, will be lost in the assembled product. However, additional nucleotides may be added between the overlap region and gene-specific sequence region to restore the pre-existing restriction site, or to introduce a new, unique restriction site. A similar principle is applied to the design of the reverse primer for the gene of interest (Fig. 4A, Right side).

Figure 4A: Assembly of Restriction Enzyme-Digested Vector and PCR-derived Insert


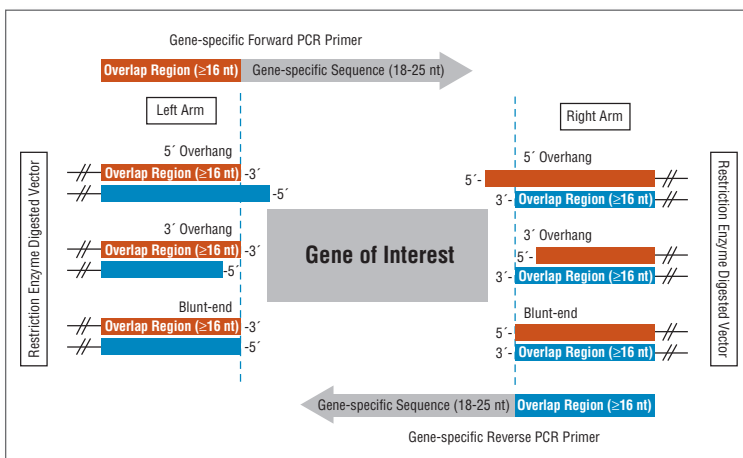


Figure 4B shows primer design for assembly of the *lacZ* gene and pMAL-c5X, digested with *NcoI* and *SbfI*. In this example, the forward primer of the gene has a "C" nucleotide (underlined) inserted between the 18-nt overlap and the N-terminal sequence of the *lacZ* gene to ensure the *lacZ* protein is in frame with the maltose binding protein.

Figure 4B: Primer Design for *lacZ* Gene and NcoI/SbfI-cut pMAL-c5X Assembly

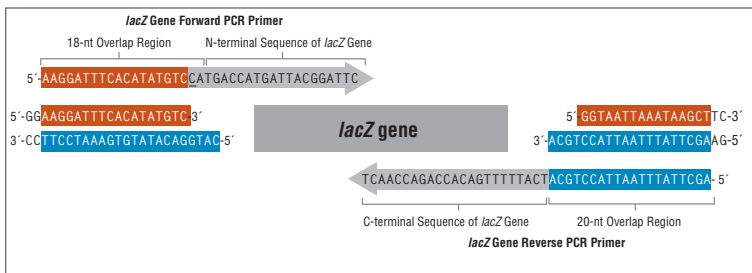

### Useful Recommendations for Vector Digestion with Restriction Enzymes

In general, the cloning vector can be linearized by any restriction endonuclease or by any combination of two restriction endonucleases displaying unique site(s) at the desired locations within the vector sequence.

*Note: Double-digestion of vector DNA with two restriction endonucleases is the best approach to reduce the uncut vector background.*

- Some restriction endonucleases cannot efficiently digest supercoiled DNA and thus may leave behind different amounts of uncut vector DNA. If not gel-purified, the uncut vector is transformable and will show up after transformation of the Gibson Assembly reaction, thereby, reducing the overall fraction of recombinant clones. A table “Cleavage of Supercoiled DNA” found at [www.neb.com/nebecomm/tech\\_reference/restriction\\_enzymes/](http://www.neb.com/nebecomm/tech_reference/restriction_enzymes/) may be used as a reference for choosing the most suitable restriction endonucleases and the number of activity units required for complete digestion of plasmid vector.
- Restriction endonucleases might have a reduced activity on plasmid DNA purified using various plasmid purification kits. In such cases, the extended restriction time or increased enzyme concentration may be necessary to digest plasmid vector to completion (or as nearly as possible to completion). When applicable, NEB highly recommends using High Fidelity (HF) restriction endonucleases to avoid star activity which may occur when digesting DNA with elevated amounts of regular restriction endonuclease for extended periods of time.
- Purification of restriction endonuclease-digested vector is not necessary unless the same restriction site is present in insert DNA. In such cases, either heat-inactivate restriction endonuclease before Gibson Assembly reaction or purify the linearized vector, either by phenol-chloroform extraction/alcohol precipitation or by electrophoresis on an agarose gel.

## Useful Recommendations for PCR Amplification

*NEB recommends using Q5 High-Fidelity DNA polymerase (NEB #M0491) or related products (NEB #M0493 or NEB #M0494) to amplify fragments of interest prior to assembly. The use of this high-fidelity DNA polymerase yields PCR products with blunt ends, thereby reducing the error rates at the fragment junctions.*

- When using circular plasmid DNA as a template, it is important to use a minimal amount of DNA (usually recommended 0.1–0.5 ng of plasmid template per 50  $\mu$ l PCR reaction) in order to reduce the template background after transformation. If higher amounts of plasmid template must be used in PCR reaction or higher amounts of PCR product must be used in the Gibson Assembly reaction, it is recommended to digest the PCR product with DpnI restriction endonuclease in order to destroy plasmid template before setting up the Gibson Assembly reaction (for protocol see below).
- Verify PCR product purity and yield by gel electrophoresis. If non-specific DNA fragments are obtained, you will need to purify the target fragment from the agarose gel to ensure the correct product assembly during the Gibson Assembly reaction.
- PCR product purification is not necessary as long as the product is > 90% pure. You can add unpurified PCR product directly from the PCR reaction into the Gibson Assembly for up to 20% of the total Gibson Assembly reaction volume (i.e. PCR products should account for 4  $\mu$ l, or less, in a 20  $\mu$ l Gibson Assembly reaction). Larger volumes of unpurified PCR products could significantly inhibit both the Gibson Assembly and the transformation. In such cases, it is recommended to column-purify PCR products and, if necessary, to concentrate DNA by ethanol precipitation.

### (Optional) DpnI Digestion Protocol

When higher amounts of plasmid template must be used in the PCR reaction, it is recommended to digest the PCR product with DpnI (NEB #R0176) restriction endonuclease in order to destroy plasmid template before setting up the Gibson Assembly reaction. DpnI cleaves only *E. coli* Dam methylase-methylated plasmid DNA, but does not cleave the PCR product since it is not methylated.

#### **DpnI Digestion Protocol:**

1. In a total 10  $\mu$ l reaction, mix 5–8  $\mu$ l of PCR product with 1  $\mu$ l of 10X NEBuffer 4 and 1  $\mu$ l (20 units) of DpnI.
2. Incubate at 37°C for 30 minutes.
3. Heat-inactivate DpnI by incubating at 80°C for 20 minutes.
4. Proceed with the Gibson Assembly Cloning procedure, described on page 11.

# Gibson Assembly Reaction:

## Optimal Quantities

NEB recommends a total of 0.02–0.5 pmols of DNA fragments when 1 or 2 fragments are being assembled into a vector. Efficiency of assembly decreases as the number or length of fragments increases. To calculate the number of pmols of each fragment for optimal assembly, based on fragment length and weight, we recommend the following formula:

$$\text{pmols} = (\text{weight in ng}) \times 1,000 / (\text{base pairs} \times 650 \text{ daltons})$$

50 ng of 5000 bp dsDNA is about 0.015 pmols.

50 ng of 500 bp dsDNA is about 0.15 pmols.

The mass of each fragment can be measured using the NanoDrop instrument, absorbance at 260 nm or estimated from agarose gel electrophoresis followed by ethidium bromide staining.

## Assembly Protocol:

1. Set up the following reaction on ice:

	Recommended Amount of Fragments Used for Assembly		
	2–3 Fragment Assembly	4–6 Fragment Assembly	Positive Control <sup>1**</sup>
Total Amount of Fragments	0.02–0.5 pmols* X $\mu\text{l}$	0.2–1 pmols* X $\mu\text{l}$	10 $\mu\text{l}$
Gibson Assembly Master Mix (2X)	10 $\mu\text{l}$	10 $\mu\text{l}$	10 $\mu\text{l}$
Deionized H <sub>2</sub> O	10-X $\mu\text{l}$	10-X $\mu\text{l}$	0
Total Volume	20 $\mu\text{l}$ <sup>***</sup>	20 $\mu\text{l}$ <sup>***</sup>	20 $\mu\text{l}$

\* Optimized cloning efficiency is 50–100 ng of vectors with 2–3 fold of excess inserts. Use 5 times more of inserts if size is less than 200 bps.

\*\* Control reagents are provided for two experiments.

\*\*\* If greater numbers of fragments are assembled, additional Gibson Assembly Master Mix may be required.

2. Incubate samples in a thermocycler at 50°C for 60 minutes. Following incubation, store samples on ice or at –20°C for subsequent transformation.

## Chemically Competent Cells Transformation Protocol:

1. Thaw chemically competent cells on ice.
2. Transfer 50  $\mu$ l of competent cells to a 1.5 ml microcentrifuge tube (if necessary).
3. If the chemically competent cells are from New England Biolabs, add 2  $\mu$ l of assembled product to NEB competent cells and go to step 4 directly. If competent cells are purchased from other manufacture, dilute assembled products 4-fold with H<sub>2</sub>O prior transformation. This can be achieved by mixing 5  $\mu$ l of assembled products with 15  $\mu$ l of H<sub>2</sub>O. Add 2  $\mu$ l of the diluted assembled product to competent cells.
4. Mix gently by pipetting up and down or flicking the tube 4–5 times. Do not vortex. Place the mixture on ice for 30 minutes. Do not mix.
5. Heat shock at 42°C for 30 seconds.\* Do not mix.
6. Transfer tubes on ice for 2 minutes.
7. Add 950  $\mu$ l of room temperature SOC media\* to tubes.
8. Place the tube at 37°C for 60 minutes. Shake vigorously (250 rpm) or rotate.
9. Warm selection plates to 37°C.
10. Spread 100  $\mu$ l of the cells onto the plates with appropriate antibiotics. Use Amp plates for positive control sample.
11. Incubate plates overnight at 37°C.

\* Please note: Follow the manufacturer's protocols for the duration and temperature of the heat shock step, as well as the optimal medium for recovery. Typically, transformation of our positive control assembly product will yield more than 100 colonies on an Amp plate with greater than 80% colonies containing inserts.

## Electrocompetent Cells Transformation Protocol:

1. Thaw electrocompetent cells on ice.
2. Transfer 50  $\mu$ l of electrocompetent cells to a pre-chilled electroporation cuvette with 1 mM gap.
3. Dilute assembled products 3-fold with H<sub>2</sub>O prior electroporation. This can be achieved by mixing 5  $\mu$ l of assembled products with 10  $\mu$ l of H<sub>2</sub>O. Add 1  $\mu$ l of the diluted assembly product to electrocompetent cells.
4. Mix gently by pipetting up.
5. Once DNA is added to the cells, electroporation can be carried out immediately. It is not necessary to incubate DNA with cells.

6. Add 950  $\mu$ l of room temperature SOC media to the cuvette immediately after electroporation.
7. Place the tube at 37°C for 60 minutes. Shake vigorously (250 rpm) or rotate.
8. Warm selection plates to 37°C.
9. Spread 100  $\mu$ l of the cells onto the plates.
10. Incubate overnight at 37°C.

## Usage Notes:

To ensure the successful assembly and subsequent transformation of assembled DNAs, NEB recommends the following:

- *Cells*: Transformation efficiency of competent cells can vary by several logs. Perceived assembly efficiency directly correlates to the competence of the cells used for transformation.
- *Electroporation*: Electroporation can increase transformation efficiency by several logs. When using the Gibson Assembly Master Mix product for electroporation, it is necessary to dilute the reaction 3-fold and use 1  $\mu$ l for transformation.
- *DNA*: Purified DNA for assembly can be dissolved in ddH<sub>2</sub>O (Milli-Q® water or equivalent is preferable), TE or other dilution buffers.
- *Insert*: When directly assembling fragments into a cloning vector, the concentration of assembly fragments should be 2–3 times higher than the concentration of vector. **For assembly of 3 or more fragments, we recommend using equimolar ratio of fragments.**
- *Biology*: Some DNA structures, including inverted and tandem repeats, are selected against by *E. coli*. Some recombinant proteins are not well tolerated by *E. coli* and can result in poor transformation or small colonies.

## Frequently Asked Questions (FAQs):

***What should I do if my assembly reaction yields no colonies, a small number of colonies, or clones with the incorrect insert size following transformation into *E. coli*?***

Assemble and transform the positive control provided with the Gibson Assembly Master Mix (see page 11,12). Successful assembly of a positive control will demonstrate that the assembly mixture is functional and the transformation conditions are suitable.

Analyze the reaction on an agarose gel. An efficient assembly reaction will show assembled products of the correct size and the disappearance of fragments.

Check the primer design of the overlapping DNA fragments to ensure that there is sufficient overlap to facilitate assembly.

Consider whether the cloned insert may be toxic to *E. coli* and a low-copy vector, such as a BAC, should be used. Because the assembled product is a covalently closed molecule, it may be alternatively amplified by PCR or RCA.

***What are the advantages of this method compared to traditional cloning methods?***

Gibson Assembly does not rely on the presence of restriction sites within a particular sequence to be synthesized or cloned. Therefore, the user has complete control over what is assembled. Furthermore, the inclusion of unwanted additional sequence, often used to facilitate the manipulation of multiple DNA sequences, can be avoided. Lastly, a greater number of DNA fragments can be joined in a single reaction with greater efficiency than conventional methods.

***How large a DNA fragment can I assemble?***

Gibson Assembly has been used to assemble and clone 300 kb DNA fragments in *E. coli*, the approximate upper limit for cloning into *E. coli*. Products as large as 1 Mb have been assembled using this approach.

***How many fragments of DNA can be assembled in one reaction?***

The number of DNA segments that can be assembled in one reaction is dependent on the length and sequence of the fragments. Gibson Assembly has been used to efficiently assemble up to twelve 0.4 kb inserts into a vector at one time. However, we recommend the assembly of five or fewer inserts into a vector in one reaction in order to produce a clone with the correct insert. A strategy involving sequential assembly can be used if all of the fragments cannot be assembled in a single reaction.



### ***How can I reduce the number of vector-only background colonies?***

To significantly reduce the background of unwanted vector-only colonies, the vector should be a PCR product rather than a restriction fragment. If background continues to be a problem, the PCR-amplified vector can be treated with DpnI to remove the template carry-over, if applicable, extracted from an agarose gel following electrophoresis.

### ***Can you PCR-amplify the assembled product?***

Yes. The assembled DNA molecule is covalently joined and may be PCR-amplified. Additionally, if the final product is a closed circular DNA molecule, it may be used as a template in rolling-circle amplification (RCA).

### ***What type of competent cells are suitable for transformation of DNA constructs created using Gibson Assembly?***

The resulting DNA constructs are compatible with most *E. coli* competent cells. NEB recommends using NEB 5-alpha Competent *E. coli* (High Efficiency, NEB #C2987). If the assembled products are larger than 10 kb, NEB recommends using NEB 10-beta Competent *E. coli* (High Efficiency, NEB #C3019) or NEB 10-beta Electrocompetent *E. coli* (NEB #C3020).

### ***Is this method applicable to the assembly of repetitive sequences?***

Yes. However, one must ensure that each DNA fragment includes a unique overlap so that the sequences may anneal and are properly assembled. The repetitive sequence can also be internalized in the first stage of a two-stage assembly strategy. If having repetitive sequences at the ends of each fragment is unavoidable, the correct DNA assembly may be produced, albeit at lower efficiency than other, unintended assemblies.

### ***What are the shortest overlaps that can be used with this assembly method?***

Productive assembly has been shown for DNA fragments with as little as 12 bp overlap. However, it depends on the GC content of the overlap. We recommend using 16 bp overlaps or more for dsDNA assembly with a  $T_m > 48^\circ\text{C}$  (AT pair =  $2^\circ\text{C}$  and GC pair =  $4^\circ\text{C}$ ).

### ***What are the longest overlaps that can be used with this method?***

The quantity of exonuclease in the Gibson Assembly Master Mix has been optimized for the assembly of DNA molecules with  $\leq 100$ -bp overlaps.

### ***Is it necessary to gel-extract restriction fragments or PCR products?***

Gel-extraction of restriction fragments is generally not necessary. A column cleanup kit or a standard phenol-chloroform extraction, followed by ethanol precipitation is sufficient. We have also used unpurified PCR products directly in assembly reactions as long as the PCR product is  $> 90\%$  pure.

***Can  $\leq 200$  bp dsDNA fragments be assembled by this method?***

Yes. For optimal results, use these fragments in  $\geq 5$ -fold excess.

***Can ssDNA oligonucleotides be combined and assembled with dsDNA fragments?***

Yes. However, the optimal concentration of each oligonucleotide should be determined. As a starting point, we recommend using 45 nM of each oligonucleotide that is less than or equal to twelve 60-base oligonucleotides containing 30-base overlaps.

***Can longer or shorter incubation times be used?***

Yes. The Gibson Assembly Kit has been optimized for assembly in 1 hour at 50°C. However, in some cases, 15 minutes may be sufficient. Reaction times less than 15 minutes are generally not recommended. Extended incubation times (up to 4 hours) have been shown to improve assembly efficiencies in some cases.

***Will the reaction work at other temperatures?***

The reaction has been optimized at 50°C, but it has been shown to work between 40°C and 50°C.

***I would like to produce overlapping dsDNA fragments by PCR. Do I need to use PCR primers that have been purified by PAGE or HPLC?***

No. Standard, desalted primers may be used.

***I would like to assemble ssDNA oligonucleotides into dsDNA fragments. Do I need to use oligonucleotides that have been purified by PAGE or HPLC?***

No. Standard, desalted primers may be used.

***Can I use a 15-nt overlap that is entirely composed of His-tag repeats (i.e. CACCACCACCACCAC)?***

No, after the His-tag, you must include at least 3 nucleotides, that are not part of the His-tag repeating sequence. Avoid repeating sequences at the end of an overlap.

***Is it necessary to inactivate restriction enzymes after vector digestion?***

No, unless the insert also carries the restriction site that was used to linearize the vector. You may still linearize the vector with such restriction enzymes, but it is necessary to heat inactivate the restriction enzyme before mixing the linearized vector with the insert in Gibson Assembly. If a heat-resistant restriction enzyme was used to linearize the vector, then vector should be purified by DNA columns, phenol-chloroform extraction or extracted from agarose gel after electrophoresis, before coming into contact with the insert.

## **Are there any differences between the Gibson Assembly Master Mix (NEB #E2611) and Gibson Assembly Master Mix included in the Gibson Assembly Cloning Kit (NEB #E5510)?**

The protocol for the Gibson Assembly Cloning Kit (NEB #E5510) is optimized to perform one to two fragment assembly and cloning into a vector, while the protocol for Gibson Assembly Master Mix (NEB #E2611) is optimized to perform multiple-fragment assemblies (up to six fragments). The major differences between the two protocols are the length of overlapping sequences between the adjacent fragments and the incubation time of the assembly reaction. The 15 minute assembly reaction protocol provided with the Gibson Assembly Cloning Kit (NEB #E5510) is recommended for assembly of fragments that are flanked by 15-25 nt overlaps. The Gibson Assembly Master Mix (NEB #E2611) is recommended for assembly of up to six fragments that are flanked by 20-80 nt overlaps, and requires a 60 minute reaction time.

## Troubleshooting

### Positive Control Yields No Colonies Following Transformation into *E. coli*

- Be sure to perform a positive control assembly reaction with 2X Gibson Master Mix, followed by transformation using competent cells from (NEB #C2987) as described on page 11,12. If competent cells are purchased from other manufacture, dilute assembled products 4 times with H<sub>2</sub>O prior transformation. This can be achieved by mixing 5 µl of assembled products with 15 µl of H<sub>2</sub>O. Add 2 µl of the diluted assembled product to competent cells.
- Competent cells may be thawed only once and cannot be repeatedly frozen and thawed without extreme loss in competency. Cells are best thawed on ice and DNA added as soon as the last bit of ice in the tube disappears.
- Do not vortex competent cells. Mix cells and DNA by gently pipetting up and down. Check cell competency by transforming 100 pg of pUC19 plasmid provided with the kit. Expect 1–3 x 10<sup>9</sup> colonies formed/µg DNA after overnight incubation on LB-ampicillin plates at 37°C.

### Gibson Assembly Reaction Yields No Colonies Following Transformation into *E. coli*

- Assemble and transform the positive control provided with the Gibson Assembly Cloning Kit. Successful assembly of a positive control will demonstrate that the Gibson Assembly Master Mix is functional and the transformation conditions are suitable.
- Check the primer design of the overlapping DNA fragments to ensure that there is sufficient and correct overlap to facilitate assembly.

- Avoid overlaps with highly palindromic sequences as they may cause up to a 10-fold reduction in recombinant colonies. When assembling fragments into a multiple cloning site (MCS) of a cloning vector, it is strongly recommended to use restriction sites that are located on the edges of the MCS to avoid overlap regions with highly palindromic sequences. Plate higher amounts of transformation reaction when using restriction sites that are located in the middle of the MCS of the cloning vector.
- Repeat the Gibson Assembly reaction using higher concentrations of fragments and/or vector. Make sure that the total volume of PCR-amplified products does not exceed 20% of Gibson Assembly reaction. If necessary, purify PCR fragments and/or PCR-amplified vector before the assembly reaction.
- Some DNA structures, including inverted and tandem repeats, are selected against by *E. coli*. Some recombinant proteins are not well tolerated by *E. coli* and can result in poor transformation.
- Test the success of the Assembly by performing PCR with primers that flank the assembled product.
- Consider whether the cloned insert may be toxic to *E. coli*, and a low-copy vector, such as a BAC, should be used. Because the assembled product is a covalently-closed molecule, it may be alternatively amplified by PCR or rolling-circle amplification (RCA).

#### Gibson Assembly Reaction Yields High Number of Clones with Incorrect Inserts

- Make sure that your PCR product is a single band of the correct size. If the PCR product is contaminated with non-specific bands, it is necessary to gel-purify the PCR product to ensure cloning of the correct insert.
- Consider whether the cloned insert may be toxic to *E. coli* and a low-copy vector, such as a BAC, should be used. Because the assembled product is a covalently closed molecule, it may be alternatively amplified by PCR or RCA.

#### Gibson Assembly Reaction Yields High Number of Small Colonies

- Some recombinant proteins are not well-tolerated by *E. coli* and can result in poor transformation efficiency or small colonies. Use a low copy number vector (i.e. pACYC184) or a vector with tight control of protein expression. When assembling into the pUC19 vector, make sure that your gene is not in frame with *lacZ* alpha fragment.

## Gibson Assembly Reaction Yields a High Number of Clones without the Insert

- PCR products may carry over large quantities of uncut plasmid template. To remove plasmid template, treat PCR products with DpnI restriction endonuclease before performing Gibson Assembly. Protocol for DpnI digestion can be found on page 10.
- Restriction enzyme digested vector may carry over large quantities of uncut plasmid. Some restriction enzymes do not cut supercoiled plasmids to completion. The best way to reduce uncut vector background is to digest the vector with two different restriction endonucleases. If a single enzyme must be used, avoid restriction enzymes that leave four-base single-stranded overhangs rich in C/G (i.e. CCGG overhang). These overhangs may self-anneal to form the transformable form of the vector molecule. Also, increase units and/or incubation time and/or purify the linear vector from agarose gel.

## References:

1. Gibson, D.G. et.al. (2009) *Nature Methods*, 343–345.
2. Gibson, D.G. et al. (2010) *Nature Methods*, 901–903.
3. Barnes, W.M. (1994) *Proc. Natl. Acad. Sci.* 91, 2216–2220.

For additional references and information on Gibson Assembly Master Mix, visit [www.sgidna.com](http://www.sgidna.com). and [www.jcvi.org](http://www.jcvi.org).

## Ordering Information

PRODUCT	NEB #	SIZE
Gibson Assembly Master Mix	E2611S/L	10/50 reactions
COMPANION PRODUCTS		
Gibson Assembly Cloning Kit	E5510S	10 reactions
Q5 High-Fidelity DNA Polymerase	M0491S/L	100/500 units
Q5 High-Fidelity 2X Master Mix	M0492S/L	100/500 rxns
Q5 Hot Start High-Fidelity DNA Polymerase	M0493S/L	100/500 units
Q5 Hot Start High-Fidelity 2X PCR Master Mix	M0494S/L	100/500 rxns
NEB 5-alpha Competent <i>E. coli</i> (High Efficiency)	C2987I/H	6 x 0.2 ml/ 20 x 0.05 ml
NEB 10-beta Competent <i>E. coli</i> (High Efficiency)	C3019I/H	6 x 0.2 ml/ 20 x 0.05 ml
NEB 10-beta Electrocompetent <i>E. coli</i>	C3020K	6 x 0.1 ml
SOC Outgrowth Medium	B9020S	4 x 25 ml





## USA

New England Biolabs, Inc.  
240 County Road  
Ipswich, MA 01938-2723  
Telephone: (978) 927-5054  
Toll Free: (USA Orders) 1-800-632-5227  
Toll Free: (USA Tech) 1-800-632-7799  
Fax: (978) 921-1350  
e-mail: [info@neb.com](mailto:info@neb.com)  
[www.neb.com](http://www.neb.com)

## CANADA

New England Biolabs, Ltd.  
Telephone: (905) 665-4632  
Toll Free: 1-800-387-1095  
Fax: (905) 665-4635  
Fax Toll Free: 1-800-563-3789  
e-mail: [info.ca@neb.com](mailto:info.ca@neb.com)  
[www.neb.ca](http://www.neb.ca)

## CHINA, PEOPLE'S REPUBLIC

New England Biolabs (Beijing), Ltd.  
Telephone: 010-82378265/82378266  
Fax: 010-82378262  
e-mail: [info@neb-china.com](mailto:info@neb-china.com)  
[www.neb-china.com](http://www.neb-china.com)

## FRANCE

New England Biolabs France  
Free Call: 0800-100-632  
Free Fax: 0800-100-610  
e-mail: [info.fr@neb.com](mailto:info.fr@neb.com)  
[www.neb-online.fr](http://www.neb-online.fr)

## GERMANY & AUSTRIA

New England Biolabs GmbH  
Telephone: +49/(0)69/305 23140  
Free Call: 0800/246 5227 (Germany)  
Free Call: 00800/246 52277 (Austria)  
Fax: +49/(0)69/305 23149  
Free Fax: 0800/246 5229 (Germany)  
e-mail: [info.de@neb.com](mailto:info.de@neb.com)  
[www.neb-online.de](http://www.neb-online.de)

## JAPAN

New England Biolabs Japan, Inc.  
Telephone: +81 (0)3 5669 6191  
Fax: +81 (0)3 5669 6192  
e-mail: [info@neb-japan.com](mailto:info@neb-japan.com)  
[www.nebj.jp](http://www.nebj.jp)

## SINGAPORE

New England Biolabs Pte. Ltd.  
Telephone: +65 6776 0903  
Fax: +65 6778 9228  
e-mail: [sales.sg@neb.com](mailto:sales.sg@neb.com)  
[www.neb.sg](http://www.neb.sg)

## UNITED KINGDOM

New England Biolabs (UK) Ltd.  
Telephone: (01462) 420616  
Call Free: 0800 318486  
Fax: (01462) 421057  
Fax Free: 0800 435682  
e-mail: [info.uk@neb.com](mailto:info.uk@neb.com)  
[www.neb.uk.com](http://www.neb.uk.com)

Some components of this product are manufactured by  
New England Biolabs, Inc. under license from  
Synthetic Genomics, Inc.

**SGIDNA**

 **NEW ENGLAND**  
**BioLabs**<sup>®</sup> Inc.  
*enabling technologies in the life sciences*