

# Synthesis and Characterization of Coordination Compounds<sup>1</sup>

Authors: D. Afzal, R. G. Baughman, H. D. Ervin, A. E. Moody, H. D. Wohlers, and J. M. McCormick\*

Previous update: March 15, 2013 & January 11, 2017; January 17, 2017 update by V. Pultz


## Introduction

Coordination compounds (also known as complex ions or simply complexes) are formed by the reaction of a Lewis acid (an electron pair acceptor, usually a transition metal) with a Lewis base (an electron pair donor), which is known as a ligand. What is unique about coordination compounds is that they are formed from chemical species that have an independent existence and that this association is often readily reversible (i. e., there is an equilibrium between the solvated metal ion and the ligand). For example,  $\text{NiCl}_2$  reacts with  $\text{NH}_3$  in aqueous solution to form the compound  $\text{Ni}(\text{NH}_3)_6\text{Cl}_2$  which contains the complex ion  $[\text{Ni}(\text{NH}_3)_6]^{2+}$ . This process is easily reversed (by the addition of  $\text{H}^+$ ) to give back the starting materials. This type of behavior was thought to be very peculiar by chemists in the 1800's. They were familiar with compounds like  $\text{CO}_2$ , which although it could be made from C and  $\text{O}_2$ , does not act like it is some loose association of C and  $\text{O}_2$ . It wasn't until the ground-breaking work of Werner (for which he won the Nobel Prize in chemistry) in the late 19<sup>th</sup> and early 20<sup>th</sup> centuries that chemists began to understand these compounds. Werner's work was greatly expanded on in the 20<sup>th</sup> century especially after it was discovered that coordination chemistry was relevant to the understanding the role of metal ions in biological systems.


You will be preparing a complex of  $\text{Co}^{3+}$  with ethylenediamine,  $\text{NH}_2\text{CH}_2\text{CH}_2\text{NH}_2$  (abbreviated: en), and a complex of  $\text{Fe}^{3+}$  with the oxalate ion,  $\text{C}_2\text{O}_4^{2-}$  (abbreviated:  $\text{ox}^{2-}$ ). Ethylenediamine and oxalate are examples of bidentate ligands, which means that they have two different atoms that can donate electron pairs to a metal ion. Ethylenediamine does this through lone pairs on its nitrogen atoms, while oxalate donates electron pairs from two of its four oxygen atoms. In these complexes the metal ion is directly bonded to six other atoms in what is called an octahedral geometry (if we connected the six atoms, the resulting solid would be an octahedron, and hence the name of this geometry). There are a number of ways in which six atoms can be arranged around a central atom in an octahedral geometry, and each of these different arrangements may give rise to compounds with the same chemical formula, but have different arrangements of their atoms (*isomers*). For example, compounds in which the actual connections between atoms (bonds) are different are called *constitutional isomers*. In this exercise you will be synthesizing and studying compounds where the bonds are the same, but the atoms are arranged differently in space (*stereoisomers*). Compounds of this type are classified as either *enantiomers* (the two compounds are mirror images of each other) or *diastereomers* (the compounds are not mirror images).

Because en is a bidentate ligand, the dichlorobis(ethylenediamine)cobalt(III) complex,  $[\text{Co}(\text{en})_2\text{Cl}_2]^+$ , that you will prepare exists as three isomers; one pair of enantiomers and their diastereomer. The isomer where the chlorides are situated on either side of the  $\text{Co}^{3+}$  ( $180^\circ$  from each other) is called the *trans* isomer (Fig. 1), while the isomer where the chlorides are next to each other in the octahedron ( $90^\circ$  from each other) is the *cis* isomer. In addition, there are two different ways in which we can put two Cl atoms *cis* to one another, and these are enantiomers


(Fig. 2). The tris(oxalato)ferrate(III) ion,  $[\text{Fe}(\text{ox})_3]^{3-}$ , exists as two enantiomers (there is no diastereomer). In one the three oxalates form a right-handed propeller, and in the other they form a left-handed propeller (both are shown in Fig. 3).

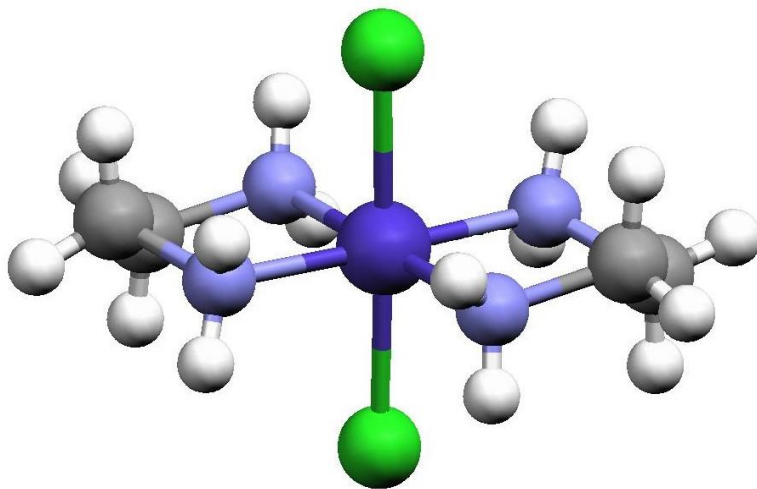

**Figure 1.** Structure of  $\text{trans}[\text{Co}(\text{en})_2\text{Cl}_2]^+$  redrawn from the [Cambridge Crystal Structure Database](#) entry CENCOS using the [Mercury](#) molecular visualization software package.

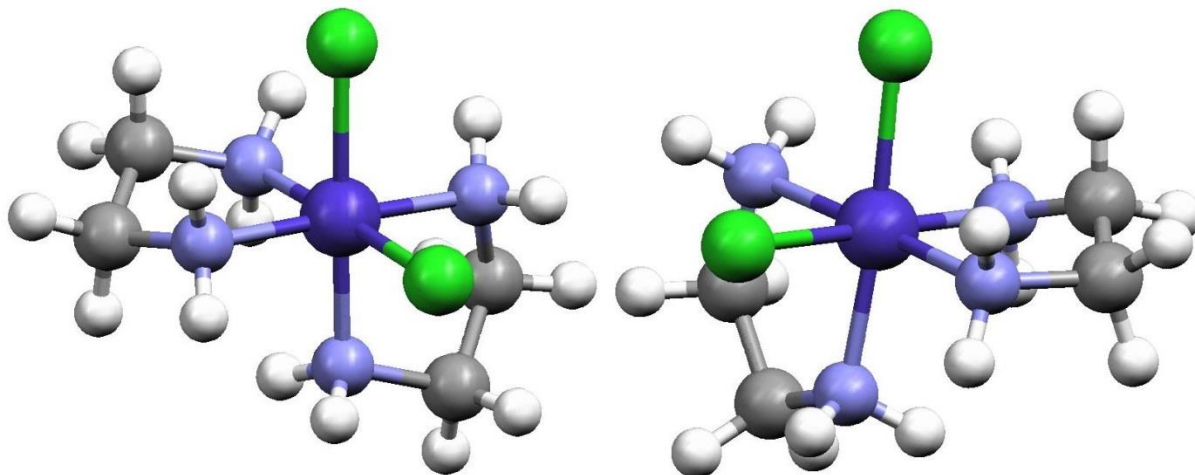

**Figure 2.** Structures of the two  $\text{cis}[\text{Co}(\text{en})_2\text{Cl}_2]^+$  enantiomers redrawn from the [Cambridge Crystal Structure Database](#) entries CENCOC and CLECO using the [Mercury](#) molecular visualization software package. The isomer on the left is designated as  $\Lambda$  (lambda) while the isomer on the right is designated as  $\Delta$  (delta) and this is determined by the orientation of the two en ligands. When one en is placed horizontally at the back of the octahedron defined by the six atoms surrounding the Co, the other en cuts across the face of the octahedron either with a positive slope ( $\Lambda$  enantiomer) or a negative slope ( $\Delta$  enantiomer).

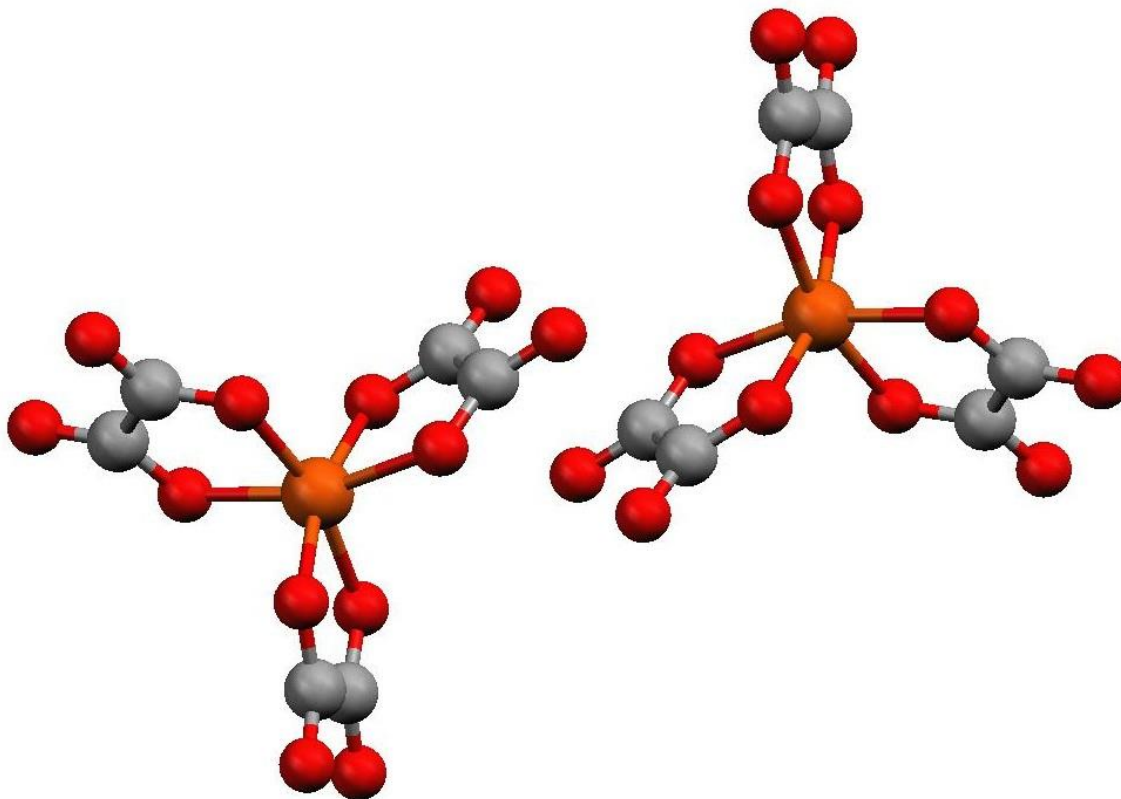


**Figure 3.** Structures of the  $[\text{Fe}(\text{ox})_3]^{3-}$  enantiomers redrawn using the [Mercury](#) molecular visualization software package from data collected by Professor Russell Baughman at Truman State University on crystals prepared by Truman students. The  $\Lambda$  enantiomer is the one on the left and the  $\Delta$  enantiomer is the one on the right. Unlike the cobalt complex, where both enantiomers can be separated and isolated by crystallization, the enantiomers of the iron complex cannot be separated. This is because there is a rapid pathway available in solution to interconvert the enantiomers.

Another interesting property of the transition metals is that their reactivity depends on several factors including the metal ion's charge, the number and type of donor atoms, and the number of d electrons present on the metal ion. It is the later property which we will exploit in this exercise both in the synthesis and then in the reactions of the compounds. The  $\text{Co}^{2+}$  ion has a  $3d^7$  electronic configuration and  $\text{Fe}^{3+}$  has a  $3d^5$  electronic configuration. Metal ions with these electronic configurations rapidly exchange their ligands and are referred to as labile. The  $3d^6$   $\text{Co}^{3+}$  is an example of an inert ion, that is, an ion that does not rapidly exchange its ligands.


In this experiment, you will synthesize  $\text{trans}-[\text{Co}(\text{en})_2\text{Cl}_2]^+$  from  $\text{CoCl}_2 \cdot 6\text{H}_2\text{O}$  according to the overall balanced chemical equation given in Scheme 1. In the mechanism by which this compound is formed  $\text{Co}^{2+}$  is initially chelated by one en and then a second en, with each step being described by an equilibrium constant. Once two, or more, en have become bound to the  $\text{Co}^{2+}$ ,  $\text{H}_2\text{O}_2$  (or, in some cases,  $\text{O}_2$  from the air) is able to oxidize the metal center to  $\text{Co}^{3+}$ . Since further substitution of a  $\text{Co}^{3+}$  complex is slow, the reaction will essentially be over at that point. Thus by careful manipulation of the amount of en present and when the oxidant is introduced, we

can force the formation of a  $\text{Co}^{3+}$  complex with either two or three en bound to the metal. Because the  $\text{Co}^{3+}$  complexes are inert we could subsequently convert  $\text{trans}-[\text{Co}(\text{en})_2\text{Cl}_2]^+$  to  $\text{cis}-[\text{Co}(\text{en})_2\text{Cl}_2]^+$  and then separate (resolve) its enantiomers.

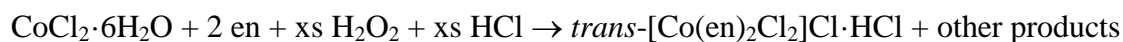


**Scheme 1.** Synthetic route to  $\text{trans}-[\text{Co}(\text{en})_2\text{Cl}_2]\text{Cl} \cdot \text{HCl}$  starting from  $\text{CoCl}_2 \cdot 6\text{H}_2\text{O}$ .

Both the *cis*- and *trans*- $[\text{Co}(\text{en})_2\text{Cl}_2]^+$  complex ions undergo hydrolysis in acidic aqueous solution, as shown in Scheme 2. This reaction can proceed through one of two mechanisms (both are shown in Scheme 3): a dissociative mechanism, where the  $\text{Cl}^-$  leaves the complex before the  $\text{H}_2\text{O}$  enters, or an associative mechanism, where the  $\text{H}_2\text{O}$  enters the complex before the  $\text{Cl}^-$  leaves.

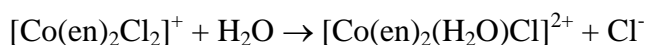


**Scheme 2.** The hydrolysis reaction of either *cis*- or *trans*- $[\text{Co}(\text{en})_2\text{Cl}_2]^+$  under acidic conditions. Note that, in general, starting with the *trans* isomer yields the corresponding *trans* isomer and starting with the *cis* isomer gives the corresponding *cis* isomer of the product.

**Associative Mechanism**

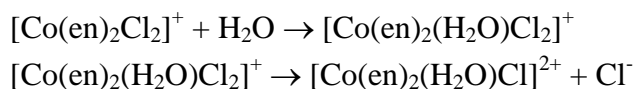


**Dissociative Mechanism**

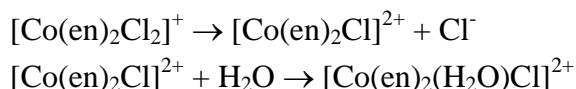

**Scheme 3.** Possible mechanisms for the acidic hydrolysis of *cis*- or *trans*- $[\text{Co}(\text{en})_2\text{Cl}_2]^+$ .


The problem is further complicated by the interconversion of the *cis*- and *trans*- isomers, as shown in Scheme 4. Luckily, this reaction is known to be much slower than the hydrolysis reaction to be studied, and as long as the reaction temperature is not taken much above  $70^\circ\text{C}$  for an extended period of time the *cis-trans* isomerization reaction need not be considered.

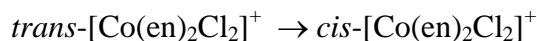

**Scheme 4.** Reaction converting  $\text{trans}-[\text{Co}(\text{en})_2\text{Cl}_2]^+$  to  $\text{cis}-[\text{Co}(\text{en})_2\text{Cl}_2]^+$ .

In the second week of this exercise, you will determine the activation energy of the first-order hydrolysis of  $\text{trans-}[\text{Co}(\text{en})_2\text{Cl}_2]^+$  under acidic conditions. You will be using a quenching method to study the reaction kinetics (compare to the direct measurement method that was used in the CHEM 130 *Kinetics of Crystal Violet Bleaching* experiment). With a quenching method a small portion of the reaction mixture is withdrawn at specific times and the reaction is then stopped by either adding reagent that stops the reaction (usually by rapidly reacting with one of the reactants) or by rapidly cooling or freezing the sample. For the reaction that we are studying, we will use the latter method to quench the reaction. Each lab group will be assigned different temperatures at which to study the reaction, and from the pooled data you prepare an Arrhenius plot and determine the reaction's activation energy.

The overall balanced chemical reaction for synthesis of the  $[\text{Fe}(\text{ox})_3]^{3-}$  is shown in Scheme 5. It is the sum of three steps, each corresponding to the stepwise reversible binding of  $\text{ox}^{2-}$  to  $\text{Fe}^{3+}$ . The metal remains in the labile 3+ oxidation state throughout the reaction, and so by adding a sufficient amount of  $\text{ox}^{2-}$ , we can force the reaction essentially completely to  $[\text{Fe}(\text{ox})_3]^{3-}$ . Because the complex is labile, we cannot resolve the  $[\text{Fe}(\text{ox})_3]^{3-}$  enantiomers, but we can use the complex's lability to confirm the 3:1  $\text{ox}:\text{Fe}^{3+}$  stoichiometry.

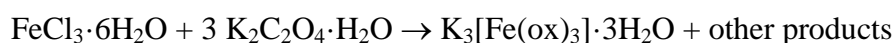

**Scheme 5.** Synthesis of  $\text{K}_3[\text{Fe}(\text{ox})_3] \cdot 3\text{H}_2\text{O}$ .

In the final week of this exercise, you will confirm the 3:1 oxalate:Fe stoichiometry in your complex in much the same way as Werner did. First you will acidify a solution of  $[\text{Fe}(\text{ox})_3]^{3-}$  which will force the sequential equilibria that formed the complex back to  $\text{Fe}^{3+}(\text{aq})$  and  $\text{H}_2\text{C}_2\text{O}_4$ . The amount of Fe and oxalate present will then be determined by redox titration against a standard  $\text{KMnO}_4$  solution.

## Experimental

For this exercise it is critical that you read each week's procedure carefully and that you and your lab partner plan what you will do before coming to the laboratory. Each week of this three-week exercise presents different challenges, and unprepared students will find it difficult to finish in the allotted time.

### Week 1

Each pair of students will need to check out the following items from the stockroom: magnetic stir bar, Büchner funnel, and filter flask.

All work of the first week's work must be carried out in the hood using the two stirring hotplates that will be in each hood. In the cobalt complex synthesis you will use one hotplate ONLY for heating, and the other ONLY for stirring. The stirring hotplate that you used for stirring will then be used in the synthesis of the iron-oxalate compound (this will allow you to start the synthesis of the iron compound while finishing the cobalt compound's synthesis). There is enough room on each plate for two beakers or two evaporating dishes; work with your hood-mates to assure that everything is done efficiently and safely.

### Synthesis of *trans*-Dichlorobis(ethylenediamine)cobalt(III) Chloride Hydrochloride<sup>2,3</sup>

Fill a 600-mL beaker approximately half full with water, and place it on a hot plate in the hood. Place your wire gauze on top of the beaker and heat to boiling.

Weigh out approximately 1 g of cobalt(II) chloride hexahydrate,  $\text{CoCl}_2 \cdot 6\text{H}_2\text{O}$ , measured to the nearest milligram. Place the cobalt salt in your evaporating dish and add 2.5 ml of distilled water. Gently swirl the dish until the cobalt has dissolved.

Once the cobalt chloride has completely dissolved in the water, add 4 mL of 10% ethylenediamine. Place your magnetic stir bar in the solution and set the evaporating dish on a cold stirring hotplate (NOT the hotplate with your water bath). Set the stir control knob so that the stir bar spins but does not splatter solution on the side of the evaporating dish (a setting of "4" on most stir plates is a good starting point). Stir for 10 min. If the evaporating dish rotates, place your triangle between the evaporating dish and the cold hotplate.

After 10 minutes, slowly and carefully add dropwise 1.6 mL of 10%  $\text{H}_2\text{O}_2$  to the solution while it is still stirring. **CAUTION!** 10%  $\text{H}_2\text{O}_2$  causes severe skin burns. Gloves are recommended and you should wash your hands after using this solution. Once the  $\text{H}_2\text{O}_2$  addition is complete, reduce the stir speed (to "3" on most of the laboratory stir plate) and stir for an additional 15 minutes.

Next, CAREFULLY add 3 mL of concentrated HCl. **CAUTION!** HCl is VERY corrosive and will cause severe burns. Immediately wash off any HCl that comes in contact with skin with copious amounts of water. Turn off the stir plate and remove the magnetic stir bar using the magnetic stir bar retriever, as your instructor will demonstrate.

Place your evaporating dish on top of the wire gauze on the beaker of boiling water. Reduce the solution's volume until there is a thick layer of dark green crystals and very little liquid left. This will take between 45 and 90 min. Monitor the level of boiling water in your water bath; you will need to gradually add water to the water bath during this time to keep it from boiling dry. While waiting, start the synthesis of potassium tris(oxalato)ferrate(III) trihydrate.

While you are waiting for the volume to be reduced, set up an ice bath. Obtain ~7 mL of methanol and put it in a small beaker or test tube in your ice bath. You will use the ice cold methanol to rinse your crystals in the final step of this synthesis.

Carefully remove the evaporating dish with the beaker tongs and set it on the bench top in the hood to cool. While your solution is cooling, set up the vacuum filtration apparatus.

Place a piece of filter paper in the Büchner funnel to collect your crystals. Carefully wet the filter paper with a couple drops of methanol BEFORE you add your crystals.

With the aspirator on, transfer the contents of your evaporating dish into the Büchner funnel using a spatula. Add about 2 mL of ice-cold methanol to your evaporating dish, swirl and pour this into the Büchner funnel. This will help remove all of your contents from the evaporating dish. Carefully pour ~5 ml of ice cold methanol over your crystals and filter. Repeat if necessary.

Carefully transfer your crystals to your cleaned evaporating dish. It is not necessary to remove the filter paper at this time. Cover with a paper towel and place it in your drawer to dry. Dispose of the liquid cobalt waste from the filtration in a designated waste container.

### **Synthesis of Potassium Tris(oxalato)ferrate(III) trihydrate<sup>4,5</sup>**

In your clean, dry 8 inch test tube dissolve 1.6 g  $\text{FeCl}_3 \cdot 6\text{H}_2\text{O}$  (measured precisely) in 4 mL of distilled water. It may be convenient to gently heat the solution in a hot water bath on a hot plate to speed the dissolution of larger chunks of the ferric chloride.

Precisely and accurately weigh out between 6.0 and 6.5 g  $\text{K}_2\text{C}_2\text{O}_4 \cdot \text{H}_2\text{O}$  and place it in a clean, dry 50-mL beaker. Add 10 mL of distilled water and gently heat on a hot plate with stirring until all of the potassium oxalate has dissolved. It is not necessary to boil the solution.

Once the oxalate salt has dissolved, quickly and carefully pour the hot oxalate-containing solution into the test tube containing the iron solution. Swirl to mix. Allow the reaction mixture to cool slowly to room temperature. Once the test tube is cool to the touch, transfer it to an ice bath and continue cooling.

Green crystals of the product may form during the initial cooling to room temperature, but they might not form until the ice cooling step. If crystals do not form after 30 min of cooling, try gently scratching the bottom of the test tube with a stirring rod or add a seed crystal (a crystal of the product obtained from another preparation), which your instructor will provide.

Decant the solution above the crystals and discard. Recrystallize the crude product by adding approximately 5 to 8 mL of hot distilled water to the crystals in the test tube, heat gently, and shake gently to affect total dissolution. Decant the green solution containing the product into a clean, dry beaker and discard any residue that remains in the test tube.

Cover the beaker with a watch glass and set it aside to cool slowly. When the beaker is cool to the touch, transfer it to an ice bath and continue to cool. Green crystals should form within about 20 min. If they do not, consult your instructor for assistance.



Collect the crystals by vacuum filtration. Wash the crystals twice with 2 mL of ice-cold distilled water and then with two 3-mL portions of acetone. **CAUTION!** Acetone is flammable, so do all these manipulations in the hood. Dry the product on the filter and then transfer it to a watch glass. Cover the watch glass with a paper towel and place it in your drawer to dry.

## Week 2

Before beginning any work this week obtain the mass of the dry *trans*-dichlorobis(ethylenediamine)cobalt(III) chloride hydrochloride and the dry potassium tris(oxalato)ferrate(III) trihydrate.

### Determination of the Activation Energy for Hydrolysis of *trans*-Dichlorobis(ethylenediamine)cobalt(III) Chloride<sup>6,7</sup>

You will use a Vernier spectrometer connected to the USB port of your laptop computer which is running LoggerPro to monitor the progress of the reaction. Operation is similar to what is described at <http://chemlab.truman.edu/instrumentation/oos/ooinstructions/> and you should quickly see the colorful display shown in Figure 2. You will make simple absorbance measurements.

At some point before beginning your kinetics runs you will need to calibrate the spectrometer using a cuvette filled with distilled water. **Important!** You must use the same cuvette for both the blank and for the actual measurements. Before inserting the cuvette into the spectrometer be sure to thoroughly wipe the cuvette's windows with a KimWipe. Any bubbles adhering to the cuvette's windows may be dislodged by **gently** tapping the cuvette with your finger. Do **NOT** tap the cuvette on the table! Be sure that the cuvette's clear windows are aligned perpendicular to the spectrometer's long side when you insert it into the spectrometer. (The spectrometers have a triangle by one side of the square hole where you insert the cuvette, and the triangle on the cuvette must also be on that side.) Also be sure that the cuvette is inserted into the spectrometer as reproducibly as possible each time. Many problems that students experience with this exercise can be traced to the cuvette not being cleaned and not being inserted properly into the instrument.

Check out a thermometer and obtain several plastic centrifuge tubes from the stockroom.

Be sure that your large test tube and all of your small test tubes are clean and dry. You will be running the reaction in the large test tube and using the small test tubes to store the aliquots removed from the reaction mixture. Place 3 ml of distilled water in each of your small test tubes and mark the water level; discard the water and dry the test tubes.

Each bench will be assigned two temperatures at which to run the reaction (the pairings are usually 40/65 °C, 45/60 °C and 50/55 °C with two benches in each lab section running the same temperature pairing, but your instructor will inform you of your actual assignments in



the laboratory). Water baths set for each of these temperatures will be in the hoods. They will be labeled and your instructor will point out which bath is in which hood. Do **NOT** change the settings on any bath! Use the thermometer to find the actual bath temperature to the nearest 0.1 °C, but do **NOT** leave the thermometer in the water bath between measurements and be careful where you set it between measurements. **Note:** it is not critical that you are exactly at your assigned temperature; the temperature reading must be stable and you must record and use the actual temperature at which the data were collected.

The actual logistics of how you and your bench mates will accomplish your assigned measurements will depend on the number of groups in the lab and the number of available spectrometers. Your instructor will give you guidance, but it is essential that you and your bench mates develop a plan of attack before coming to the laboratory.

Place 30 mL of 0.01 M HNO<sub>3</sub> in your large test tube and place it in the water bath. Place 1 mL of cold distilled water in a small plastic centrifuge tube (volumes are marked on the side). Keep the centrifuge tube cold in an ice bath until ready for use.

To perform a kinetics run, precisely weigh out approximately 0.2 g of *trans*-[Co(en)<sub>2</sub>Cl<sub>2</sub>]Cl·HCl. Place this in the small centrifuge tube containing the cold water, close the tube and shake vigorously to dissolve the solid. Pour the entire contents of the centrifuge tube into your large test tube and mix thoroughly. Do this as quickly as possible; it is best if this can be accomplished without removing the large test tube from the water bath. Start timing from when the solution is completely mixed.

Remove 3 mL of the reaction mixture immediately upon mixing. Place it in one of your small test tubes and place the tube in the ice bath. Obtain the absorbance spectrum of this sample after it has cooled back to about room temperature and you have time. Record the absorbance at 505 nm and at 800 nm (this is your data point at  $t = 0$ ). Don't forget to thoroughly wipe the cuvette's windows with a KimWipe before you place it in the spectrometer. Any bubbles adhering to the cuvette's windows may be dislodged by **gently** tapping the cuvette with your finger. Do **NOT** tap the cuvette on the table! If an absorbance measurement at 800 nm is not approximately zero, about 0.020 or smaller, re-wipe the cuvette's windows and try the measurement again.

Repeat this process until you have removed a total of nine samples, but make sure ~3 mL remains for the 70-°C water bath. The timing between samples will vary with temperature. At 40 °C you must follow the reaction for at least 40 minutes, but at 65 °C you will only need to follow the reaction for 5 minutes. Work out the timing of the sample withdrawals before you start a run. **You must know precisely how long each sample was in the water bath after complete mixing until it was placed in a small test tube which was immediately put in an ice bath.** You will use times and absorbances for a graph.


Once you have removed nine samples, place the test tube containing the remaining 3 mL of the reaction mixture in the 70-°C water bath for no more than five minutes. Remove the test tube from the bath and place it in the ice bath. Once it has cooled to near room temperature, obtain the reaction mixture's absorbance spectrum, and record the absorbance at 505 nm and 800 nm; this is your  $t = \infty$  data point.

Prepare a fresh sample and obtain data at the other assigned temperature. Repeat the measurements at the two temperatures as time permits.

All cobalt-containing waste is to be placed in the proper container. Any unused cobalt complex will be saved for later use. Please put it in the container provided, as directed by your instructor.

### *Week 3*

Use the method of half-reactions to balance the two redox reactions that you will use in this analysis **before** coming to the lab.

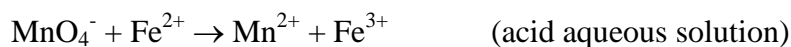

The optimal amount of titrant is around 20.00 mL. In this experiment the titrant has an approximate molarity of 0.01 M  $\text{MnO}_4^-$  and you use very pure  $\text{K}_2\text{C}_2\text{O}_4 \cdot \text{H}_2\text{O}$  to determine the molarity more precisely. Based on the optimal amount of titrant and its approximate molarity, how much pure  $\text{K}_2\text{C}_2\text{O}_4 \cdot \text{H}_2\text{O}$  should be your target mass? Show all work.

Prepare titration tables using

[http://vpultz.sites.truman.edu/files/2014/06/Hydrolysis-of-Oil-of-Wintergreen\\_titration-tables.pdf](http://vpultz.sites.truman.edu/files/2014/06/Hydrolysis-of-Oil-of-Wintergreen_titration-tables.pdf)

as a model, but make appropriate changes.

### **Standardization of 0.01 M Potassium Permanganate Solution**

Obtain approximately 75 mL of un-standardized 0.01 M  $\text{KMnO}_4$  solution. **CAUTION!**

Potassium permanganate is a strong oxidizing agent! Immediately wash any spilled on your skin with copious amounts of water.

Wash your buret several times, as your instructor will demonstrate, with no more than 2 mL of the permanganate solution. Each time allow the solution to drain through the buret's tip into a waste beaker. After the final wash fill the buret to near the 0.00-mL mark such that the upper part of the meniscus is at or below the 0.00-mL mark. Make sure that there are no bubbles in the buret tip. Run some of the solution through to clear any bubbles (gentle taping might help dislodge any recalcitrant bubbles). Refill the buret, if necessary. Record the initial volume of solution in the buret. Because the  $\text{MnO}_4^-$  ion is so strongly colored, make all your volume readings from the top of the meniscus, and your instructor may allow you to record to the nearest 0.1 mL.

Accurately weigh out the approximate mass of  $\text{K}_2\text{C}_2\text{O}_4 \cdot \text{H}_2\text{O}$  that you calculated you will need to standardize the  $\text{MnO}_4^-$  solution (helpful hint: it should be around 0.1 g) and record the actual mass. Place the oxalate in an Erlenmeyer flask and add about 30 mL of distilled water and 5 mL of 6 M  $\text{H}_2\text{SO}_4$ . Swirl the flask gently to dissolve the solid.

Place the flask on a stirring hotplate in the hood and carefully heat the solution to approximately 60 °C. Do **not** boil! Remove the flask from the hotplate (a folded paper towel makes an excellent hot pad). This titration must be performed fairly quickly so that the solution does not cool too much. Be sure that you swirl the flask as you add the permanganate solution and wash down any solution that splashes on the side of the flask with a small stream of distilled water. If the solution cools too much, reheat. Titrate the sample with the permanganate solution until the slightest purple color persists for at least 30 sec. Record the final buret reading. Calculate the volume of permanganate used and from it determine the  $[\text{MnO}_4^-]$ . Repeat until three determinations agree with each other within 2%.

### **Determination of Oxalate in Potassium Tris(oxalato)ferrate(III) Trihydrate**

Accurately weigh out about 0.1 g of your  $\text{K}_3[\text{Fe}(\text{ox})_3]\cdot 3\text{H}_2\text{O}$  and record the mass to the nearest 0.001 g. Put in an Erlenmeyer flask. Add 30 mL distilled water and 5 mL of 6 M  $\text{H}_2\text{SO}_4$ . Swirl to dissolve the solid and then heat to 60 °C, again taking care not to boil the solution. Titrate as described above to the first permanent purple color. Record the buret readings, calculate the volume of titrant dispensed, and determine the %  $\text{C}_2\text{O}_4^{2-}$  by mass in the sample.

Do **NOT** discard the solution at this point; proceed to the iron analysis given below.

### **Determination of Iron in Potassium Tris(oxalato)ferrate(III) Trihydrate**

Take your faintly-purple sample from the oxalate analysis above, and set in on a hotplate in the hood (there may be a small amount of brown precipitate forming at this point, but do not be concerned). Gently heat it until it almost boils and add approximately 100 mg of Zn to the hot solution. Cover the flask with a watch glass and continue heating until the yellow color (from  $\text{Fe}^{3+}$ ) disappears ( $\text{Fe}^{2+}$  is colorless in solution). While you are waiting, do more oxalate titrations and set up a gravity filtration with fluted filter paper (ask your instructor to demonstrate how to flute filter paper, if you are unsure).

Quickly filter the hot, colorless solution into another Erlenmeyer flask, again using a folded paper towel to handle the hot flask. It is important that this filtration be done quickly to minimize the amount of  $\text{Fe}^{2+}$  that is re-oxidized to  $\text{Fe}^{3+}$  by  $\text{O}_2$  in the air. Rinse the funnel with several small (approximately 5 mL total volume) portions of distilled water. Titrate this solution with the permanganate solution in your buret until the first faint trace of a persistent pink color. Record initial and final buret readings. Calculate the volume of permanganate used and determine the % Fe by mass in your sample.

Repeat the two titrations until you run out of sample, or when three of the oxalate titrations agree with each other within 2%.

## Results and Analysis

### Week 1

Show your calculations for the theoretical yield of *trans*-[Co(NH<sub>2</sub>CH<sub>2</sub>CH<sub>2</sub>NH<sub>2</sub>)<sub>2</sub>Cl<sub>2</sub>]Cl·HCl and K<sub>3</sub>[Fe(C<sub>2</sub>O<sub>4</sub>)<sub>3</sub>]·3H<sub>2</sub>O. Assume that CoCl<sub>2</sub>·6H<sub>2</sub>O and FeCl<sub>3</sub>·6H<sub>2</sub>O are the limiting reagents. After you have determined the mass of your dry *trans*-[Co(en)<sub>2</sub>Cl<sub>2</sub>]Cl·HCl and K<sub>3</sub>[Fe(ox)<sub>3</sub>]·3H<sub>2</sub>O, calculate a percent yield for each compound. Show your calculations.

### Week 2

[Prepare a graph in Excel](#) showing absorbance as a function of wavelength in which you overlaid all of the spectra for one run. The scale on the y-axis should be 0 to 1 (you may need to set the maximum value on the y-axis to 1.1 or 1.2, depending on your actual concentrations) and the x axis scale should be set from 400 to 800 nm. If the data were obtained correctly, the spectra should all have an absorbance of about 0 at 800 nm and there should be two *isosbestic* points (points where the absorbance does not change as a function of time, indicating that only two species were contributing to the absorbance throughout the reaction).

From the absorbance at 505 nm at each time subtract the absorbance at 800 nm at the same time; this corrects for any baseline drift or other variations that occurred between samples. With the corrected absorbance readings at 505 nm ( $A_t$ ) and your corrected absorbance at infinite time ( $A_\infty$ ), prepare a first-order integrated rate law graph (i. e.,  $\ln(A_\infty - A_t)$  as a function of time) for each run. We use  $A_\infty - A_t$  instead of concentration in these graphs because 1) we are following the formation of the product over time and 2) the reactant has a relatively large absorbance at this wavelength.

From each graph determine the value of the rate constant at that temperature. Report each rate constant (along with its uncertainty [at 95% confidence](#)) and the temperature at which it was obtained to the instructor and the rest of the class. From the class data (rate constant as a function of temperature), prepare an Arrhenius plot, determine the activation energy and its confidence interval at the 95% confidence limit.

Does the fact that the hydrolysis reaction is first-order in the cobalt complex and your value of the activation energy allow you to conclusively rule out one of the possible mechanisms? You will need to derive the rate law predicted by each mechanism for each of the possible cases: 1) the first step is irreversible and rate determining, 2) the first step rapidly goes to equilibrium and the second step is slow, 3) the first step is reversible, but does not come to equilibrium and the second step is slow. What additional experiments could you do to help distinguish between the possible mechanisms?

### Week 3

Show all raw data and calculated values in titration tables. Include sample calculations for trial 1 in each of the three tables. Start with standardization of permanganate. Calculate the average percent oxalate by mass in your K<sub>3</sub>[Fe(ox)<sub>3</sub>]·3H<sub>2</sub>O and share this with the class. From the class data calculate the average percent oxalate by mass in K<sub>3</sub>[Fe(ox)<sub>3</sub>]·3H<sub>2</sub>O, the [standard deviation](#) and the [confidence interval at the 95% confidence limit](#). Be sure to perform a [Q-test](#) on the class data first to exclude any suspect point. Calculate the true percent oxalate by mass and determine a percent error for both your data and the class data.

Calculate the average percent iron by mass in your  $\text{K}_3[\text{Fe}(\text{ox})_3] \cdot 3\text{H}_2\text{O}$ . Share this with the class, and calculate the average percent iron by mass in  $\text{K}_3[\text{Fe}(\text{ox})_3] \cdot 3\text{H}_2\text{O}$ . Perform a Q-test on the class data first to exclude any suspect point. Calculate the true percent iron by mass and determine a percent error for both your data and the class data.

## Conclusions

This exercise contains elements of both a [synthesis](#) exercise and a measurement exercise

## Summary of Results

In your *Summary of Results* you should have tables that summarize the results for each week: Week 1, the percent yield of *trans*- $[\text{Co}(\text{en})_2\text{Cl}_2]\text{Cl} \cdot \text{HCl}$  and  $\text{K}_3[\text{Fe}(\text{ox})_3] \cdot 3\text{H}_2\text{O}$ ; Week 2, the average rate constants at each temperature that you determined and the activation energy that was determined from the class data (along with a 95% confidence interval for each, if possible); and Week 3, the statistical (average, standard deviation and confidence interval) results for the  $\text{K}_3[\text{Fe}(\text{ox})_3] \cdot 3\text{H}_2\text{O}$  analyses.

## References

1. <http://chemlab.truman.edu/chemical-principles/>
2. Szafran, Z.; Pike, R. M. and Singh, M. M. *Microscale Inorganic Chemistry: a Comprehensive Laboratory Experience*; John Wiley and Sons: New York, 1991.
3. Bailar, J. C., Jr. *Inorg. Synth.* **1946**, 2, 222-225.
4. Beran, J. A. *Laboratory Manual for Principles of General Chemistry, 5<sup>th</sup> Ed.*; John Wiley and Sons: New York, 1994.
5. Marcus, S.; Sienko, M. J. and Plane, R. A. *Experimental General Chemistry*; McGraw-Hill: New York, 1988.
6. Jolly, W. L. *Encounters in Experimental Chemistry, 2<sup>nd</sup> Ed.*; Harcourt, Brace, Jovanovich: New York, 1985, 77-79.
7. Herrick, R. S.; Mills, K. V. and Nestor, L. P. *J. Chem. Educ.* **2008**, 85, 1120-1122. Click [here](#) to view a PDF version of this article (Truman addresses and *J. Chem. Educ.* subscribers only).